

MP Biomedicals, LLC

FastPrep-24™ 5G Instruction Manual (v2.7)



[This page intentionally left blank]

Copyright

All rights reserved. No part of this book may be reproduced, stored in a retrieval system, or transmitted, in any form or by any means (e.g., electronic, mechanical, photocopying, recording, or otherwise) without the prior written permission of MP Biomedicals, LLC.

Disclaimer

Information in this document is subject to change without notice. MP Biomedicals, LLC, makes no guarantees of any kind with regard to this material, including but not limited to; the implied warranties or merchantability and fitness for a particular purpose. MP Biomedicals, LLC shall not be held liable for errors contained herein or for incidental or consequential damage in connection with the furnishing, performance, or use of this material.

Manufacturer

MP Biomedicals, LLC
Corporate Headquarters
3 Hutton Centre Drive, Suite 100
Irvine, CA 92707, USA

MP Biomedicals United States
29525 Fountain Parkway
Solon, OH 44130, USA

Customer Service / Technical Support

Tel: 800-854-0530

Fax: 800-334-6999

Email: custserv@mpbio.com, biotech@mpbio.com

Website: www.mpbio.com

Part Number: 116005500

Revision 2

Rev Date, September 14, 2017

Document No.: TEC-FP24-002

1	SECTION I: INTRODUCTION	7
2	A. General Description	7
3	B. Principle of Operation	10
4	C. FastPrep-24™ 5G Features	11
5	D. FastPrep-24™ 5G Specifications	12
6	SECTION II: SAFETY	14
7	A. Regulatory	14
8	B. Safety	14
9	C. Environmental Conditions	15
10	D. Biological Safety	15
11	SECTION III: INSTALLATION	16
12	A. Unpacking	16
13	1) To unpack the FastPrep-24™ 5G:	17
14	2) Inspection	17
15	3) Additional Materials Required:	18
16	B. Installation	19
17	1) Connecting the Power	20
18	2) Set-up, Controls and Functions	22
19	3) Installation Verification	24
20	SECTION IV: OPERATION	25
21	A. Powering up the FastPrep-24™ 5G	25
22	B. Dome Operation	25
23	C. Loading and Securing the Samples, Adaptor and Cam Lock Installation	26
24	D. Operating the FastPrep-24™ 5G	39
25	E. 'Emergency Stop Button' Operation	39
26	SECTION V: SOFTWARE	40
27	A. FastPrep-24™ 5G 'Touch Screen' Software Operation	40

TABLE OF CONTENTS

28	1) Running a Recommended Program Assay:	40
29	2) Manually Programming an Assay:	44
30	F. Running an Assay	46
31	G. Saving an Assay	48
32	3) Using Saved Programs:	49
33	A. Retrieving a Saved Program	49
34	B. Deleting a User-Programmed Assay:	50
35	4) Stopping Cycle in Progress:	51
36	B. FastPrep-24™ 5G Administrative Functions	52
37	1) About FastPrep-24™ 5G: System Versions and End-User Agreement	53
38	2) Date /Time Set	54
39	3) Optional Settings	55
40	A. Keyboard Clicks	56
41	B. Buzzer Warnings	56
42	C. Strobe Effect	57
43	4) Software Update	57
44	a) Installing the FastPrep-24 5G Host Application	58
45	b) Install FastPrep-24 5G driver	58
46	c) Open the FastPrep-24 5G in Safe Mode	61
47	d) Install the Firmware Update	62
48	e) Calibrating the FastPrep-24™ 5G 'Touch Screen	65
49	5) Exporting History File	65
50	6) Technical Support	68
51	<i>Resetting the FastPrep-24™ 5G:</i>	69
52	SECTION V: WARRANTY & LIABILITY	70
53	SECTION VI: MAINTENANCE	71
54	A. Maintenance	71
55	B. Yearly Maintenance	Error! Bookmark not defined.
56	C. Fuse Replacement (as needed)	73
57	SECTION VII: PARTS INFORMATION	75
58	APPENDIX 1: FASTPREP® ADAPTORS	83
59	LARGE-SCALE ADAPTORS	83

TABLE OF CONTENTS

60	HIGH THROUGHPUT ADAPTORS	84
61	CRYOGENIC ADAPTORS	85
62	METAL ADAPTORS	86
63	APPENDIX 2: FASTPREP® LYSING MATRIX	87
64	APPENDIX 3: RECOMMENDED PROGRAM SETTINGS	89
65	APPENDIX 4: FASTPREP® INSTRUMENTS FOR OTHER APPLICATIONS	94
66	FastPrep-96™ Instrument	94
67	SuperFastPrep-2™ Portable Homogenizer, 2 x 2 mL, battery operated	96
68		

SECTION I: INTRODUCTION

A. General Description

The FastPrep-24™ 5G (FP5G) Instrument is a high-speed, benchtop reciprocating instrument for the optimal lysis of challenging and routine sample types. It is intended for use in applications that require grinding, lysing or homogenization of various solid sample materials. Sample types include but are not limited to the following: human, animal and plant tissues; bacterial, yeast and fungi cells teeth, bones; and soil, fecal samples as well as mineral or building material.

The FastPrep-24™ 5G is a software controlled, standalone instrument, designed with a user-friendly touch screen interface as well as unique features such as an increased speed and a strobe light to indicate adapter rotation. The FastPrep-24™ 5G comes standard with the QuickPrep-3 adapter, which allows for the simultaneous processing of 24 samples in 2 mL tubes; additional adapters are available for different size tubes.

Samples are placed into tubes along with a specific, optimized lysing matrix, which are subsequently placed into the appropriate adapter within the instrument. The FastPrep-24™ 5G will partially rotate the adapter, and therefore the tubes, in a unique manner, resulting in a multidirectional, simultaneous beating of the lysing matrix on the sample material. The beating action will result in a mechanical lysing of the cellular membrane, allowing for the extraction of the target material (i.e., DNA, RNA, or protein) into the solution. The speed and time of the agitation, in addition to the adapter used, will vary based on specific needs, and individual parameters can be programmed using the FastPrep-24™ 5G touch screen.

The touch screen visual display allows for user-friendly control of the instrument. The FastPrep-24™ 5G, while programmed with predefined assays designed for the optimal lysis of routinely run sample types, also allows for the programming and saving of custom assays.

A front and back view of the FastPrep-24™ 5G Instrument, along with each of the major components, is illustrated in [Figure I-1](#) and [Figure I-2](#) respectively.



Figure I-1: Major Components of the FastPrep-24™ 5G (Front View)

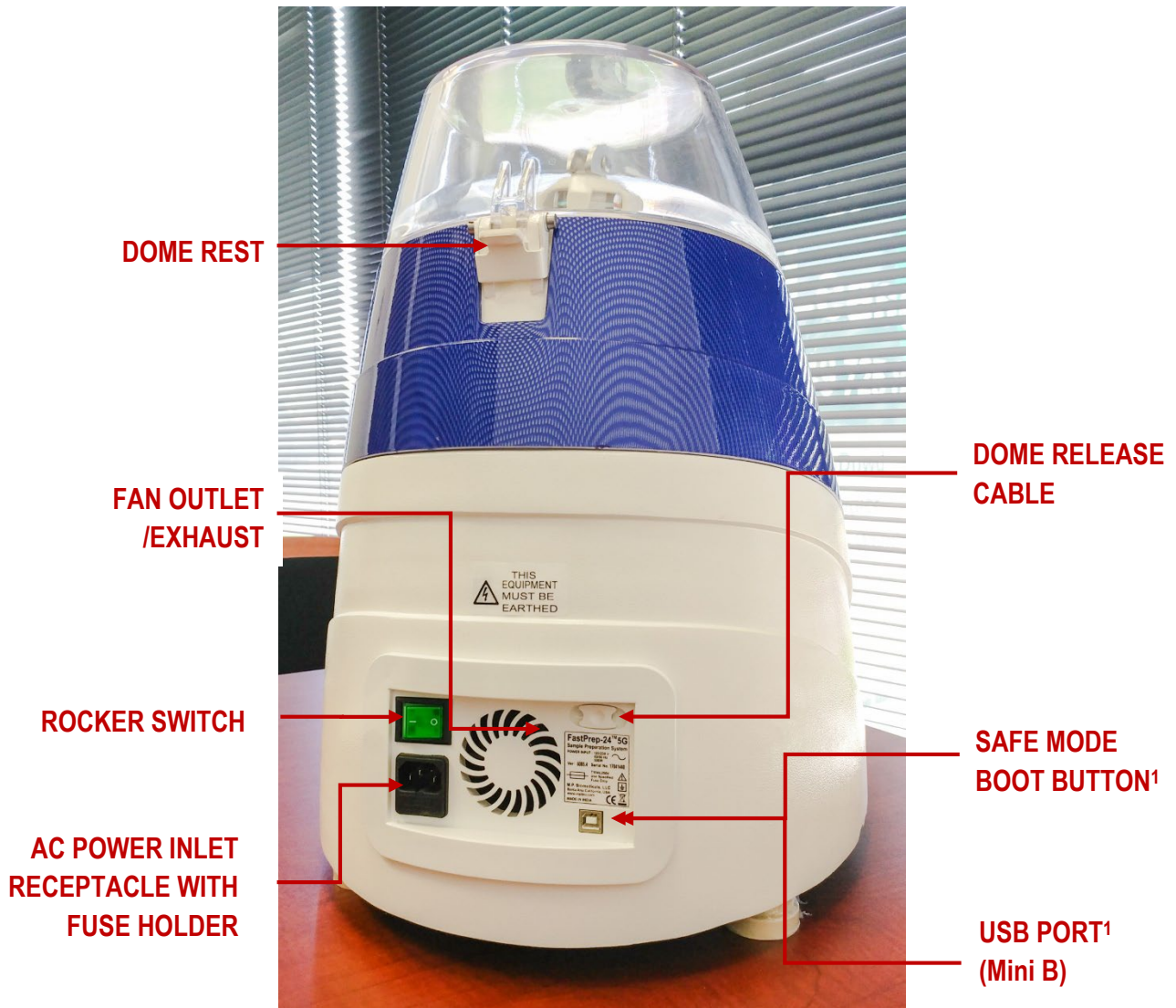


Figure I-2: Major Components of the FastPrep-24™ 5G (Back Panel View)

¹ Required for software update

B. Principle of Operation

The FastPrep-24™ 5G is a high-speed, benchtop instrument intended for use in a sample preparation application where it is necessary to grind, lyse, or homogenize difficult and routine samples, including but not limited to, biological tissues, fluids and excretions; cell cultures; environmental samples; and other inorganic solid matrices.

The FastPrep-24™ 5G causes cellular disruption in samples by facilitating the collision of the sample with a sample specific lysing matrix, resulting in the mechanical disruption of the cellular membrane and release of the target analyte into the surrounding protective buffer. The FastPrep-24™ 5G is designed to maximize the quality and quantity of the yield through efficient cellular disruption, and the effective removal of inhibitors and removal or inhibition of nucleases and proteases.

The FastPrep-24™ 5G facilitates efficient and effective cellular lysis using Lysing Matrix tubes and a reciprocating motion. Individual samples are placed into individual tubes along with the appropriate lysing matrix; the tubes are sealed and disposable, minimizing the potential for contamination. The tubes are placed directly into the samples holder or adapter housed within the instrument, which produces a reciprocating motion that results in the collision of the sample with the lysing matrix. The rate of collision and energy of impact, critical factors in cellular disruption efficiency, are a function of the FastPrep-24™ 5G instrument speed setting and the specific gravity of the lysing matrix. The rate of collision is proportional to the speed, while the energy of impact is proportional to the square of the speed. The FastPrep-24™ 5G offers variable speed and time settings for optimal cellular disruption from a wide variety of sample types.

C. FastPrep-24™ 5G Features

The FastPrep-24™ 5G offers a variety of optional, interchangeable Adaptors to hold tubes of various sizes, allowing for versatility in processing volumes as well as lysis under cryogenic conditions. The FastPrep-24™ 5G comes standard with the QuickPrep-3 adaptor, allowing for a maximum load of 24, 2 mL samples, processed at ambient temperatures. Additional Adaptors available, as well as the maximum load and volume for each, are indicated below. Detailed information on each adaptor can be found in [Appendix 1](#).

Description	Catalog No.	Sample Load	Weight Empty (g)	Weight Filled (g)
QuickPrep-3	116005512	24 x 2 mL	175	264
BigPrep	116002525	2 x 50 mL	570	749
TeenPrep	116002526	12 x 15 mL	506	806
HiPrep	116002527	48 x 2 mL	467	643
CoolPrep (Cryogenic)	116002528	24 x 2 mL	680	768
CoolTeenPrep (Cryogenic)	116002530	6 x 15 mL	627	757
CoolBigPrep (Cryogenic)	116002531	2 x 50 mL	700	815
TallPrep	116002540	24 x 4.5 mL	417	596
All-Metal QuickPrep	116002545	24 x 2 mL	307	396
All-Metal TeenPrep	116002546	12 x 15 mL	840	1140
All-Metal BigPrep	116002547	2 x 50 mL	500	684



DAMAGE TO INSTRUMENT

Use of Adaptors other than those listed in this operations manual may damage the FastPrep-24™ 5G

D. FastPrep-24™ 5G Specifications

The FastPrep-24™ 5G has the following specifications:

FastPrep-5G Instrument Specifications	
Item	Specification
Product Name:	FastPrep-24™ 5G
Model:	6005
Software:	Microcontroller: ARM Processor Memory: 64 MB App Version (Touch Screen): 1.3.2 Firmware Version (Driver): 1.3.7.2 Programming: Field Alterable Flash, 12 End-User Programmable Assays
Interface:	Touch Screen Interface Screen Size: 151.5 mm x 85.5 mm
Controls:	Programmable Settings for Speed, Adapter, Time, Lysing Matrix, Sample Quantity, Sample Units, Cycle, Pause Time; Display Readout
Programmable Assays:	12 Programmable Assay Locations
Pre-Defined Assays:	72 Pre-defined and Optimized Assay Programs
Time Range:	1 to 120 Seconds in 1 Second Increments (Default: 40 Seconds)
Speed Range:	4 to 10 m/sec in 0.5 m/sec Increments; (Default: 6 m/sec for QuickPrep-3 Adapter)
Cycles:	1 to 9 Cycles (Default: 1 Cycle)
Adapter(s):	Standard QuickPrep-3, 24 x 2 mL Adaptor; Additional Adaptors Available
Pause Time:	1 to 300 Second Pause Between Cycle in 1 Second Increments (Default: 300 Seconds)
Lysing Matrix:	Lysing Matrix A through Z, Plus Others
Quantity:	Free Entry
E-Stop:	Switch Push DPST-NC (10A, 110V)
Date / Time Set:	Available as 24-Hour Clock

FastPrep-5G Instrument Specifications	
Data Export:	Via USB
Acceleration:	< 2 seconds to maximum speed
Deceleration:	< 2 second to stop
Dimensions:	Height: 490 mm Base: 472 mm x 385 mm (Elliptic Shape)
Weight:	23.6 kg (52 lb)
Power Requirement:	120 VAC / 60 Hz, 6 A 230 VAC / 50 Hz, 3 A
Protecting Fuse:	T4A 205V, 5x30 mm
Operating Air Temperature:	2 to 48 °C (35 to 118 °F)
Operating Relative Humidity:	30 to 55%
Overvoltage Category:	II
Maximum Sound Level:	< 70 dB
Maximum Altitude:	2000 m (6562 feet)

SECTION II: SAFETY




A. Regulatory

The FastPrep-24™ 5G is US Class 1 device. It is compiling with USFDA general control and CE marked with the essential requirements of the applicable European laws or Directives with respect to safety, health, environment, and consumer protection.

B. Safety

The FastPrep-24™ 5G is designed for safe operation. However, for your safety and the safety of others, the FastPrep-24™ 5G should be operated as specified in this Operations Manual. Failure to do so could result in injury to yourself or others, or in damage to the instrument. There are various symbols on the FastPrep-24™ 5G unit relating to safety, which are indicated in the table below.

Safety Symbols used with the FastPrep-24™ 5G

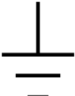Symbol	Description
	Alternating Current
	Resistor
	Electrical Equipment, Dispose Properly
	Caution, Warning
	Power On
	Power Off
	Earth Terminal Ground

C. Environmental Conditions

The FastPrep-24™ 5G is designed to be safe under the following environmental conditions:

FastPrep-24™ 5G Operating Specifications

Environmental Condition	Specification
Location:	Indoor use only
Maximum altitude:	2000 m
Maximum sound level:	70 dB
Overvoltage Category:	II
Temperature:	2 to 48 °C (35-118°F)
Humidity:	< 95% RH
Ventilation	≥ 5 cm around instrument
Voltage	110 VAC/60 Hz or 230 VAC/50H Hz

D. Biological Safety

The FastPrep-24™ 5G may be used with samples that contain materials of human origin. All blood or other potentially infectious materials should be considered infectious regardless of the perceived status of the source individual. Observe Standard Precautions when working with potentially infectious materials. Ensure all materials are disposed of in accordance with applicable regulations.



DANGER

Some samples used with the FastPrep-24™ 5G may contain materials of human origin. Observe Standard Precautions when working with potentially infectious materials.

SECTION III: INSTALLATION

A. Unpacking

MP Biomedicals will ship the FastPrep-24™ 5G directly to your location. Inspect the exterior shipping crate for damage upon arrival and notify MP Biomedicals immediately in case of damage.

The FastPrep-24™ 5G shipping box has been labeled to indicate the following:



PRECISION INSTRUMENT

Handle instrument with care.



Keep package upright.

Store the FastPrep-24™ 5G shipping crate upright in a secure location until ready to install. For installation purposes, the FastPrep-24™ 5G shipping crate is oriented as indicated on the box.



HEAVY OBJECT

The FastPrep-24™ 5G should be lifted with both hands firmly under the base of the instrument, using the finger-locator cut-outs provided.



The FastPrep-24™ 5G must be transported in its original packaging. Use of packaging other than the original packaging may damage the instrument and/or void the product warranty.

INSTALLATION**1) To unpack the FastPrep-24™ 5G:**

1. Open top flaps of shipping container.
2. Remove the smaller pieces of protective foam and air-bag stuffing from inside the box. The FastPrep-24™ 5G will be packaged within two large foam pieces.
3. Firmly grip each piece of foam, one with each hand, and lift foam and instrument vertically from the box. The foam pieces are designed to slide easily out of the box while cradling the instrument safely. Alternatively, lay the box on its side and grip the foam pieces and slide the instrument out.
4. Place the foam and instrument combination on a sturdy table or bench top.
5. Remove each piece of foam from around the instrument, freeing the instrument.
6. Place all packaging materials back into the box and store for future use.

NOTE: Alternately, the FastPrep-24 outer container may be placed horizontally on a bench top or floor for foam and instrument removal.

2) Inspection

The items listed in the [“Packing List”](#) below are included with the FastPrep-24™ 5G as provided by MP Biomedicals, LLC. Please contact MP Biomedicals, LLC., if an item is not received as listed.

FastPrep-24™ 5G Packing List

Part Description	Quantity
FastPrep-24™ 5G	1 ea
Power cord (European)	1 ea
Power cord (US)	1 ea
USB Cable	1 ea
Replacement Spring/Screw Assembly	2 ea
Cam Lock	1 ea

Fuse 10 Amp	2 ea
Strike Bolt	2 ea
Replacement QuickPrep-3Adaptor	1 ea
FastPrep-24™ 5G Operations Manual	1 ea

3) **Additional Materials Required:**

The FastPrep-24™ 5G, when shipped to the end user facility, contains all necessary materials for immediate instrument use following proper installation, as well as some backup materials such as fuses. However, during the lifetime of the product, items such as fuses may require replacement more often than the backups provided allow for. Use of the FastPrep-24™ 5G with parts other than those specified by the manufacturer may impact performance.

- 5 x 20 mm Miniature Fuse, Time-Lag T, L, 250 VAC

The FastPrep-24™ 5G contains the different components as indicated in [Figure I-1](#) and [Figure I-2\(Section I\)](#). Inspect the FastPrep-24™ 5G unit for damage that may have been caused by shipping. In the event of damage or questions, please contact MP Biomedicals, LLC.

B. Installation

The FastPrep-24™ 5G should be installed in an indoor location only, within an operating environment that meets all specifications as indicated. The FastPrep-24™ 5G should be placed on a clean, dry, stable surface within 1.2 m of an electrical outlet, and in a location that allows for a minimum of 5 cm space around the instrument base at all times to ensure proper ventilation.

1. Place the FastPrep-24™ 5G on a sturdy table or bench top.
2. Position the FastPrep-24™ 5G in its final installation position, ensuring easy access to the power cord and receptacle.
3. Disengage the 'Emergency Stop' by turning clockwise until it pops out. **(Device is shipped with 'Emergency Stop' engaged for safety and will not run until disengaged.)** ([Figure III-1](#))
4. To open the 'Dome', press ANY of the four options on the FastPrep-24™ 5G 'Home Screen' ([Figure V-1](#)) and press [OPEN]. Alternatively, pull the 'Dome Release Cable' on back plate of device to release the 'Dome Latch' ([Figure III-2](#)). Raise the 'Dome' gently until it reaches the rest position. For more on 'Dome' operation see ['Dome' Operation](#).
5. Remove any remaining foam from the chamber assembly.
6. Gently lower the 'Dome' to a closed position. Engage the 'Dome Latch' by pressing downward on the 'Dome' until the latch mechanism clicks.



Figure III-1: Disengage 'Emergency Stop'



Figure III-2: 'Manual Dome Release Cable'

1) Connecting the Power

The power cord connects to the back of the FastPrep-24™ 5G as seen in [Figure III-3](#). The FastPrep-24™ 5G can operate on either 100-120 VAC/60 Hz or 220-240 VAC/50 Hz power. Ensure that the main 'Rocker Switch' in the back of the instrument is in the OFF position. Connect the power cord to the instrument power fixture (below 'Rocker Switch') and then connect the other end to a compatible power supply outlet.



Figure III-3: Power Cord Connection

To ensure a secure connection, the FastPrep-24™ 5G should be located no more than 1.2 m from an appropriate electrical outlet.

NOTE:

Ensure the 'Rocker Switch' (O/ I), located on the rear panel, is in the OFF [O] position prior to connecting the power supply.



IMPORTANT

For safe use, ensure that the FastPrep-24™ 5G is properly grounded by connecting the power cord to an approved electrical outlet.

2) Set-up, Controls and Functions



IMPORTANT

To ensure safe operation and optimal performance of the FastPrep-24™ 5G instrument, read this Operations Manual before operating the instrument.

The FastPrep-24™ 5G instrument comes fully assembled and requires very little set-up once installed. Set-up of the FastPrep-24™ 5G, and control of the instrument during operation, is accomplished through the use of a 'Touch Screen' based central control panel as indicated in [Figure III-4](#).



Figure III-4: FastPrep-24™ 5G Touch Screen Control Panel

Users interface with the FastPrep-24™ 5G using a 'Touch Screen', which contains the different menu functions. The [main menu functions](#) of the FastPrep-24™ 5G are indicated below.

FastPrep-24™ 5G Touch Screen Main Menu Functions

Item Name	Operation	Function
Display screen	<i>Display</i>	Displays menus and programmed settings; counts down run times
Recommended Programs	<i>Press</i>	Allows user to access pre-defined program by sample type (i.e., plants, bacteria, environmental, human/animal, yeast & fungi, and feces)
Saved Programs	<i>Press</i>	Allows user to access saved custom programs
Program Manually	<i>Press</i>	Allows user to program custom parameters (i.e., speed, adapter, time, lysing matrix, quantity and unit, cycles and pause time)
Admin Tools	<i>Press</i>	Allows user to access administrative functions such as date/time set, export data, options, etc.

The FastPrep-24™ 5G has [custom programmable settings](#) as indicated:

Variable Speed and Time Settings

Item	Range	Increment
Speed	4.0 to 10.0 ¹ m/s	0.5 m/s
Adapter	All FastPrep-24 Adapters	See adapter list (Appendix 1)
Time	0 to 120 ² sec	1 sec
Lysing Matrix	All MP Lysing Matrix Types	See Appendix 2
Quantity	Free entry	NA
Quantity Unit	mg, g, mL, cells	NA
Cycles	1 to 9	1
Pause Time	0 to 300 sec	1 sec

NOTE:

Speed settings higher than 6.5 m/s are available only for the QuickPrep adapters. All other optional adapters are limited to a maximum speed of 6.5 m/s.

IMPORTANT

While the FastPrep-24™ 5G is capable of run times up to 120 seconds, **it is highly recommended that any run time over 60 seconds be monitored very closely!** Some Lysing Matrix combinations can generate significant heat after 60 seconds at high speeds and tube failure is possible.

3) Installation Verification

Prior to using the FastPrep-24™ 5G for the first time, an installation verification should be performed. To verify installation:

1. Load and secure the QuickPrep-3 'Adaptor' as described in [Section IV, OPERATION.](#)
2. Run the FastPrep-24™ 5G at 6 m/sec rpm for 60 seconds (Information on programming the FastPrep-24 may be found in [Section V, SOFTWARE](#)). Listen for loud noises, grinding or whining noises.

IMPORTANT

If loud noises, grinding or whining noises can be heard with the FastPrep-24™ 5G, immediately engage the ['EMERGENCY STOP BUTTON'](#). Contact MP Biomedicals, LLC.

3. If the 6 m/sec run is acceptable, run the FastPrep-24™ 5G at 8 m/sec for 60 seconds. Listen for loud noises, grinding or whining noises.
4. If the 8 m/sec run is acceptable, run the FastPrep-24™ 5G at 10 m/sec for 60 seconds. Listen for loud noises, grinding or whining noises.
5. Turn instrument off when not in use by using the 'Rocker Switch' located on the back panel.

NOTE:

See Section V for [Administrative Functions](#) such as setting [\[DATE/TIME\]](#), [Optional Settings](#), and [End-User Agreement and Registration](#) of your new FastPrep-24™ 5G Instrument.

SECTION IV: OPERATION

A. Powering up the FastPrep-24™ 5G

Turn on the FastPrep-24™ 5G by switching the 'Rocker Switch' to the **[ON]** position ([Figure I-2](#)). The FastPrep-24™ 5G 'Touch Screen' will light up, and the instrument will display 2 splash screens of MP Biomedicals information, followed by the FastPrep-24™ 5G 'Home Screen' ([Figure V-1](#)).

NOTE:

Ensure that the 'Emergency Stop Button' has been pulled all the way out. The unit will not start if the 'Emergency Stop Button' is engaged ([Figure III-1](#)).

**IMPORTANT**

The samples and 'Adaptor' must be secured properly before the instrument is operated. Failure to secure either could result in personal injury or damage to the instrument.

B. Dome Operation

The safety 'Dome' is locked and unlocked by an electromechanical latch assembly (Dome latch), which is operated via the 'Touch Screen' user interface.

1. To open the Dome, press any of the four options on the FastPrep-24™ 5G 'Home Screen'.
2. Press [OPEN]. The mechanism will activate and release the 'Dome Latch'.
3. Carefully lift the 'Dome' and slowly rotate backwards on its hinge until it contacts the 'Dome Rest'.
4. To close the 'Dome', carefully lower the 'Dome' back into the closed position and press firmly downward at the 'Dome Latch' until the 'Dome Latch' automatically engages.

NOTE:

The [OPEN] button is available on all programming screens, but is not available on some Admin screens or when a run is in progress.

NOTE:

An assay may be programmed but cannot be initiated with the 'Dome' open.

NOTE: In the event of power loss, or to open the ‘Dome’ manually, pull the ‘Dome Release Cable’ on the back panel of the device ([Figure III-2](#)).

C. Loading and Securing the Samples, Adaptor and Cam Lock Installation

The FastPrep-24™ 5G comes standard with a QuickPrep-3 ‘Adaptor’. The maximum sample load for a FastPrep-24™ 5G, using the QuickPrep-3 ‘Adaptor’, is 24 each 2 mL tubes. Alternative ‘Adaptors’ will have different sample load specifications; information on the alternative ‘Adaptors’ available is provided in [Appendix 1](#) and information on the Lysing Matrices used with each ‘Adaptor’ can be found in **Appendix 1**. For optimal performance, MP Biomedicals recommends the use of FastPrep® Lysing Matrix with the FastPrep-24™ 5G instrument.



IMPORTANT

Some ‘Adaptors’ must be appropriately balanced for safe operation.



IMPORTANT

Use Standard Precautions when handling potentially infectious samples.


For convenience, MP Biomedicals offers many different Lysing Matrices for both challenging and routine sample types.

NOTE: *It is strongly recommended that only Lysing Matrix Tubes from MP Bio be used with the FastPrep-24™ 5G instrument.* While tubes from other suppliers may be compatible, MP Bio Lysing Matrix Tubes are designed with optimal dimensions and strength. Improper tube dimensions may result in lower lysing efficiency, increased chance of tube failure, and potential damage to ‘Adaptors.’

Once the samples are prepared, the samples are loaded onto the QuickPrep-3 Adaptor. The QuickPrep-3 Adaptor is then loaded onto the FastPrep-24™ 5G as follows:

1. At the FastPrep-24™ 5G ‘Home Screen’ ([Figure V-1](#)), press any of the 4 options.

2. Press the [OPEN] button to release the 'Dome Latch'.
3. Lift the 'Dome' and allow to rest in the open position.
4. Loosen the 'Cam Lock' by lifting handle upward and around the 'Handle Pivot', until the 'Cam Lock' is in the unlocked position ([Figure IV-1.1](#)). This requires moderate force and it may be necessary to steady the 'Adaptor' with free hand. ([Figure IV-1](#)).
5. Rotate the 'Cam Lock', in the unlocked position, counterclockwise to loosen until the threads disengage.

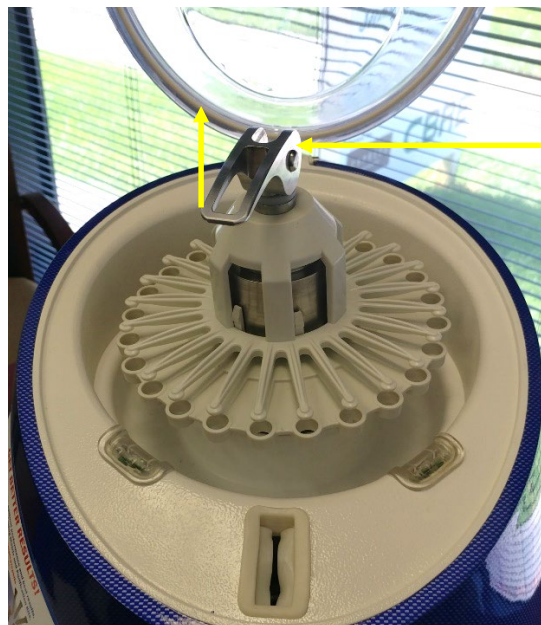

**CAM LOCK IN
'LOCKED' POSITION**

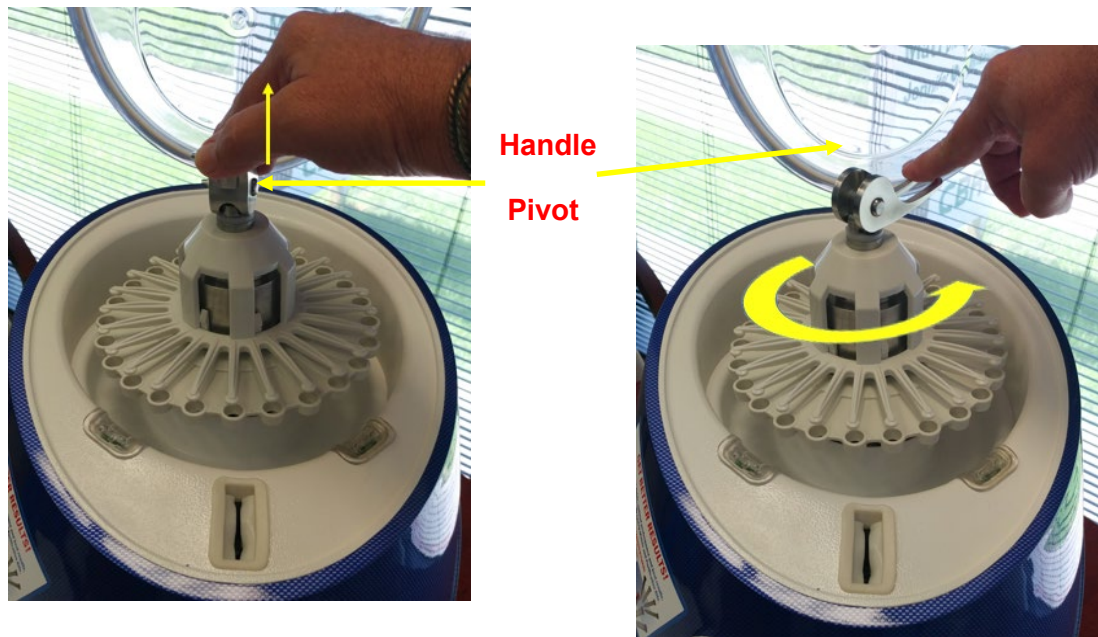


Figure IV-1: QuickPrep-3 and Unlocked Cam

6. Remove the 'Cam Lock' from device.
7. Remove the 'Adaptor' assembly ([Figure IV-2](#)).

CAM LOCK LOCKED POSITION




CAM LOCK UNLOCKED POSITION



Figure IV-1.1: Cam Lock Locked or Unlocked



Spoke Plate

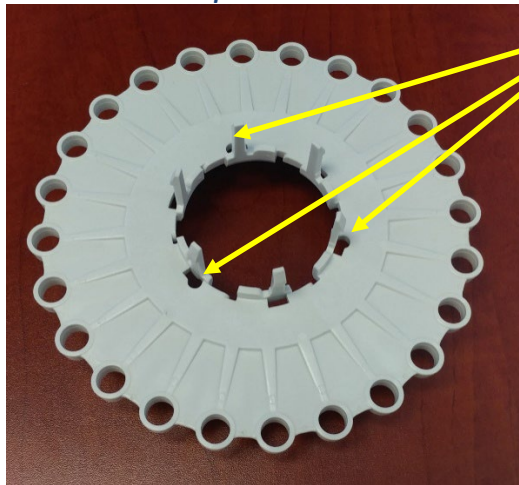

**THREE (3)
ALIGNMENT HOLES
THAT ACCEPT
LOCKING PIN (SEE
STEP #11)**

Tube Deck



Assembled, Unlocked Position

Figure IV-2: QuickPrep-3 Assembly Components

- Lift the 'Spoke Plate' slightly and rotate clockwise so that the retention spokes move away from the holes on the 'Tube Deck', leaving them open for loading ([Figure IV-3](#)).



Figure IV-3: QuickPrep-3 Sample Loading Position

- Load the sample tubes into the holes of the 'Tube Deck', and press so that they seat snugly and the 'Tube Base-Ring' contacts the 'Tube Deck' surface ([Figure IV-3.1](#)). For less than full loads, tubes must be balanced symmetrically.

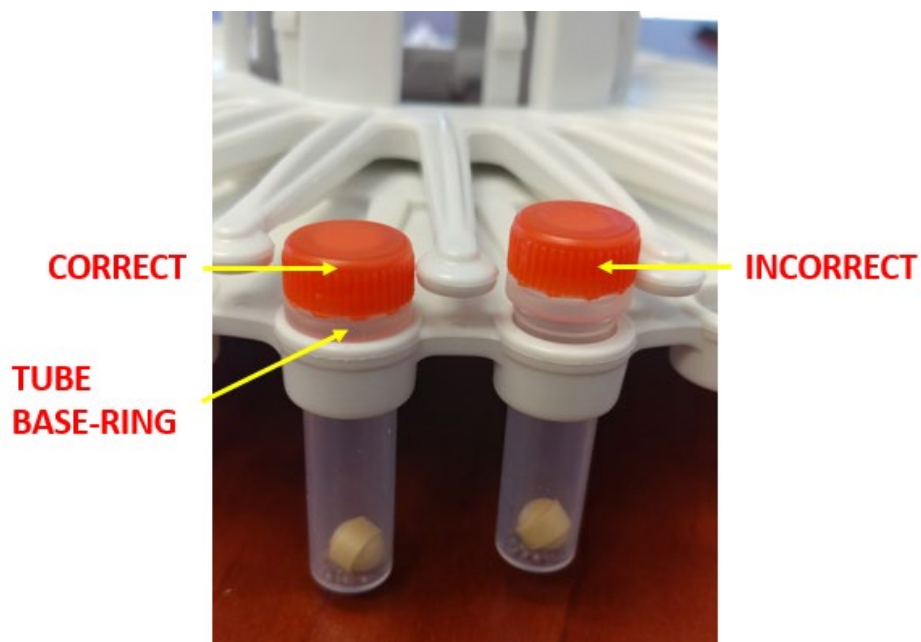


Figure IV-3.1: Tube Loading

10. Lift the 'Spoke Plate' slightly again and rotate it counter-clockwise so that the retention spokes are placed directly above each sample tube cap ([Figure IV-4](#)).



Figure IV-4: Spoke Plate Placement

11. When all the sample tubes have been loaded, place the 'Adaptor' back into the FastPrep-24™ 5G. Align the locking pin of the 'Three Step Assembly' with an alignment hole under the 'Adaptor', ensuring its proper placement ([Figure IV-5](#)). **Rotate the 'Adaptor' on the '3-Step Assembly' until the locking pin engages, no more than $\frac{1}{3}$ turn, and the adaptor will drop down on the pin ~5 mm ([Figure IV-6](#)).**



WARNING

Failure to properly position and secure the 'Adaptor' within the FastPrep-24™ 5G may cause damage to the 'Adaptor' and/or the instrument.

NOTE:

Refer to ([Figure IV-2](#)) and ([Figure IV-5](#)) for views of 'Alignment Holes' 'Three Step Assembly' and 'Locking Pin.'

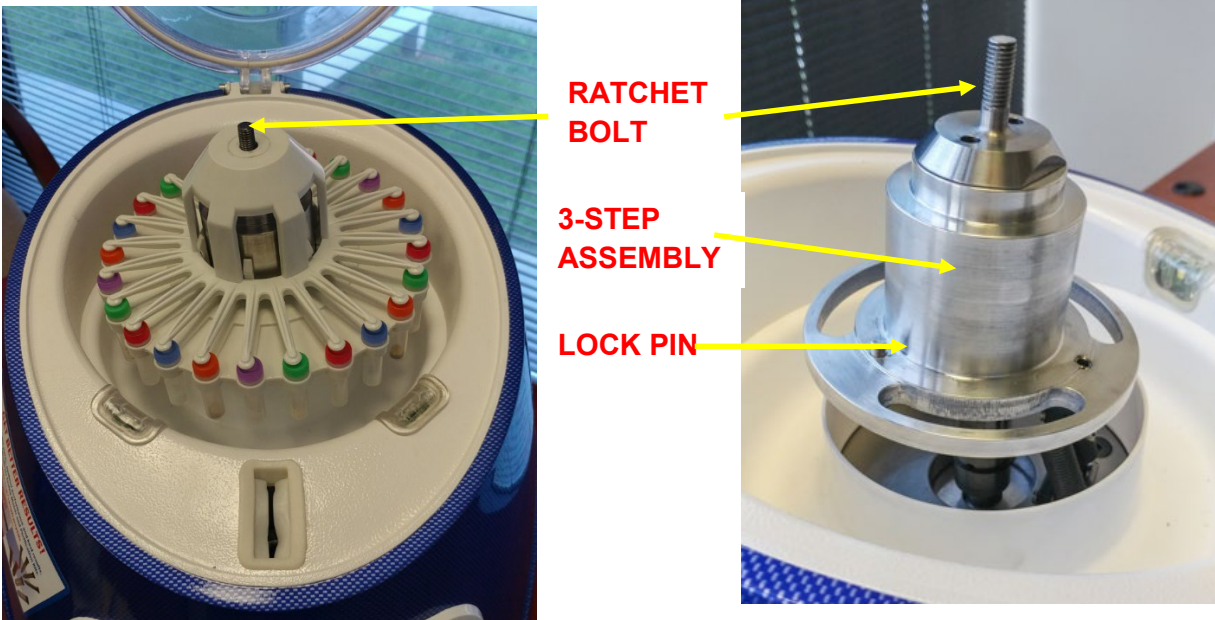


Figure IV-5: Adaptor Placement

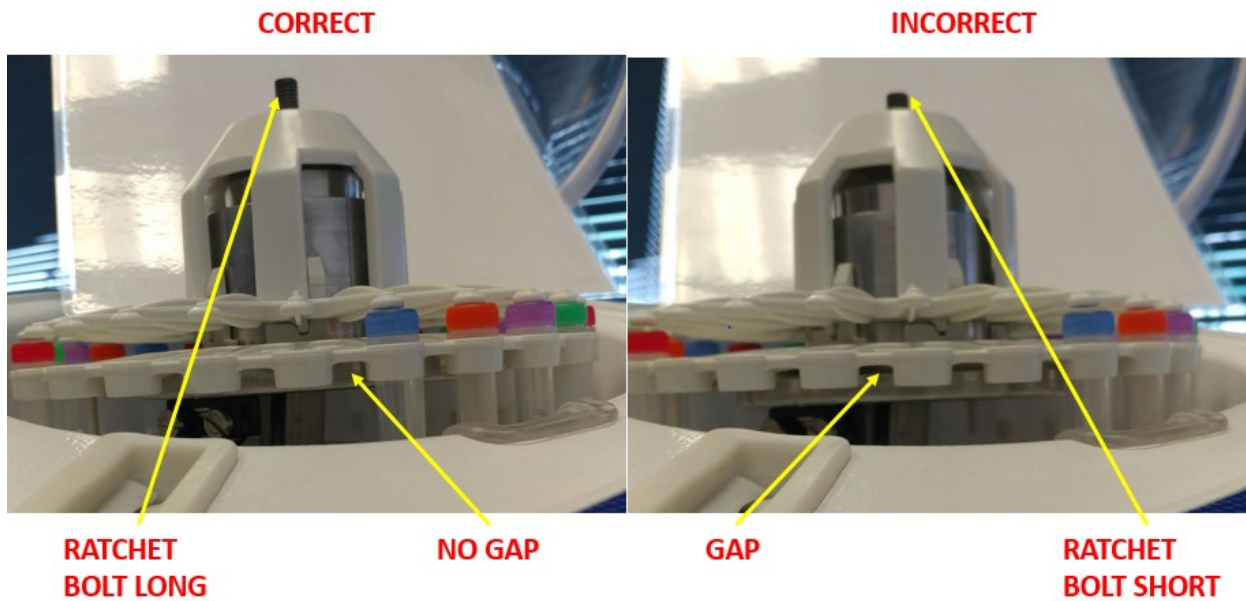


Figure IV-6: Rotate Adaptor until Seated

12. Place the 'Cam Lock' in the locked position onto the threaded 'Ratchet Bolt' and turn clockwise to engage and seat threads ([Figure IV-7](#)).

**CAM LOCK
LOCKED
POSITION**



Figure IV-7: Thread Cam Lock on Ratchet Bolt

13. Rotate 'Cam Lock' clockwise to tighten. There should be no resistance and rotation should be done using one finger ([Figure IV-8](#)).

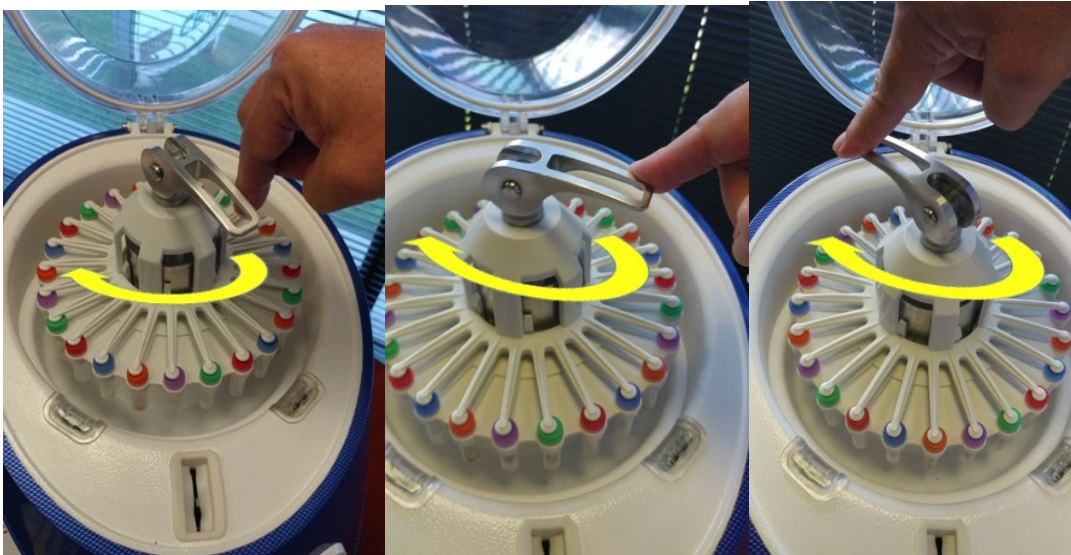


Figure IV-8: Rotate Cam Lock Clockwise

14. After ~ 3 complete clockwise revolutions, the 'Cam Lock' base will contact the 'Adaptor' and resistance will be felt. Stop tightening at this point.

15. Lift the 'Cam Lock' handle straight up and around the 'Handle Pivot' to achieve the unlocked position. [\(Figure IV-9\)](#) Make sure that the 'Cam Lock' **DOES NOT ROTATE** around the 'Ratchet Bolt' while unlocking.

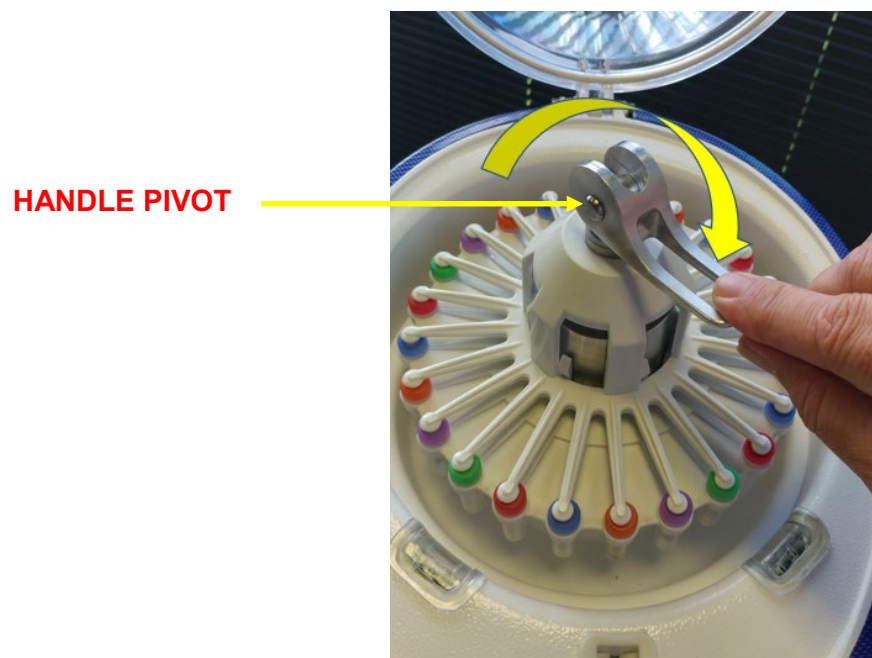


Figure IV-9: Unlock Cam Lock, 180 Degrees about Pivot

16. Rotate the unlocked 'Cam Lock' ~ 180 degrees (~ ½ turn) clockwise around the 'Ratchet Bolt' [\(Figure IV-10\)](#). The degrees' rotation may vary slightly and should be verified and adjusted, if necessary, See Step #17.



Figure IV-10: Rotate Unlocked Cam Lock 180 Degrees

17. Keep the 'Cam Lock' in the same position with respect to the 'Ratchet Bolt' and lift the handle upwards and around the 'Handle Pivot' to lock. Use free hand to steady the 'Adaptor.' [\(Figure IV-11\)](#). As the handle is lifted, no resistance will be felt at first. **Resistance will be felt when the handle reaches vertical, or slightly past vertical. If resistance is felt before vertical, loosen the 'Cam Lock' slightly by turning a few degrees counterclockwise about the 'Ratchet Bolt.'** **Never turn 'Cam Lock' on 'Ratchet Bolt' in the locked position with resistance. Always unlock it before adjustments are made.**



WARNING

Tightening or loosening the 'Cam Lock' on the 'Ratchet Bolt' in the Locked Position with Resistance (Pressure) may cause damage to the 'Cam Lock' and/or the instrument.

NOTE:

It is recommended to steady the 'Adaptor' with your free hand to allow a stronger grip to lock the 'Cam Lock' as well as to ensure no rotation around the 'Ratchet Bolt' occurs. See [\(Figure IV-11\)](#) and [\(Figure IV-12\)](#).

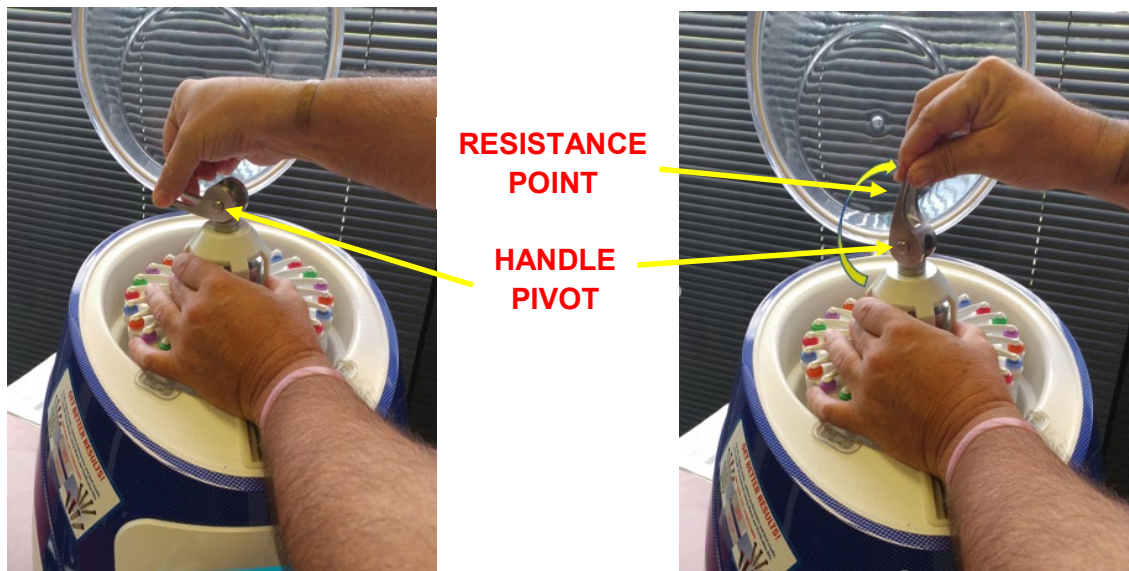


Figure IV-11: Lift Handle Around Pivot

18. Continue to pull the handle over the 'Handle Pivot' and press it downward, with force, until it stops. [\(Figure IV-12\)](#). It may be necessary to re-grip the handle.



Figure IV-12: Press Handle Down Until Stop

19. "Cam Lock" is now fully locked. [\(Figure IV-13\)](#) Tubes and 'Adaptor' are properly positioned and secure.



Figure IV-13: Cam Lock Fully Locked

20. Close the 'Dome' and press down firmly until the 'Dome Latch' engages.

NOTE:

The FastPrep-24™ 5G 'Dome' contains an electromechanical lock mechanism. The 'Dome Latch' must be engaged for the instrument to operate.

WARNING



When first installing the 'Cam Lock' system, carefully observe the instrument motion once the assay is started. If properly installed, 'Cam Lock' will remain fixed in place while the motor runs. If any movement of the 'Cam Lock' is observed (loosening, counterclockwise rotation) immediately engage the [\(EMERGENCY STOP BUTTON\)](#) and repeat the 'Cam Lock' installation procedure. Failure to stop the motor could result in personal injury or damage to the instrument

21. Once sample processing is complete, to remove 'Adaptor' from instrument, first unlock 'Cam Lock' by lifting handle upward and around the handle pivot ([Figure IV-14](#)), until the 'Cam Lock' is in the unlocked position ([Figure IV-15](#)). This requires moderate force. Steady the 'Adaptor' with free hand for better grip and to eliminate any rotation around 'Ratchet Bolt.'

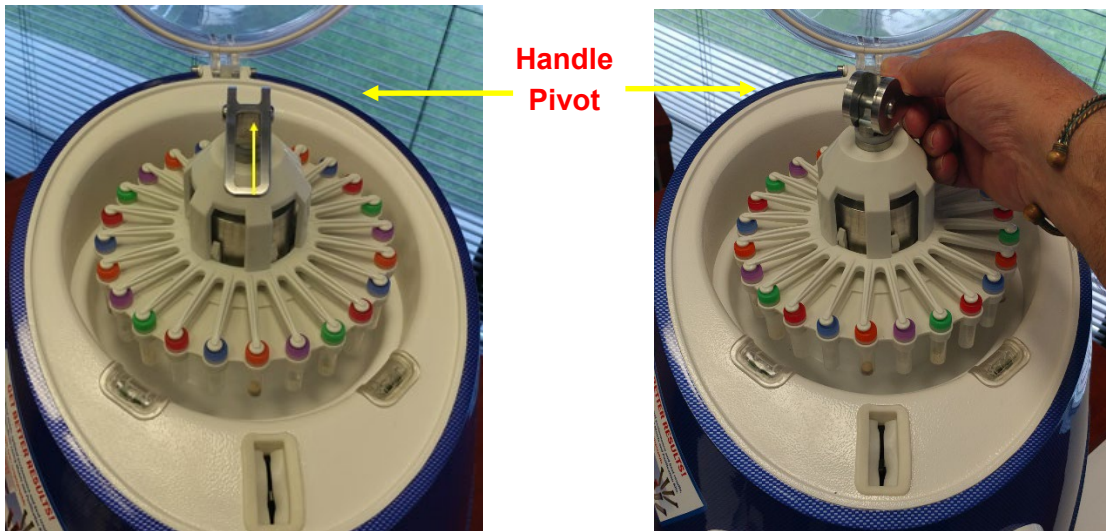


Figure IV-14: Lift Cam Handle Upwards Figure IV-15 Cam Handle Unlocked

22. Rotate the 'Cam Lock', in the unlocked position, counter-clockwise to loosen until the threads disengage ([Figure IV-16](#)). Remove 'Cam Lock' from device ([Figure IV-5](#)).



Figure IV-16 Unscrew Unlocked Cam Lock

23. 'Adaptor' can now be removed from device.

D. Operating the FastPrep-24™ 5G

The FastPrep-24™ 5G is primarily operated using the FastPrep-24™ 5G custom software, accessed through the 'Touch Screen' interface. For details on the FastPrep-24™ 5G software operation, refer to [Section V: SOFTWARE](#).



IMPORTANT

The samples and 'Adaptor' must be secured properly before the instrument is operated. Failure to do so could result in personal injury or damage to the instrument.

E. 'Emergency Stop Button' Operation

The red 'Emergency Stop Button' ([Figure IV-17](#)), located directly below the 'Touch Screen' is a safety feature and can be used at any time. Engaging the 'Emergency Stop Button' will immediately cut all power directed to the FastPrep-24™ 5G instrument.

1. To engage the 'Emergency Stop Button', press the button firmly until it clicks. This will cut all power to the FastPrep-24™ 5G.
2. To restore power, firmly rotate the 'Emergency Stop Button' clockwise until it releases.



Figure IV-17: 'Emergency Stop Button'

SECTION V: SOFTWARE

A. FastPrep-24™ 5G 'Touch Screen' Software Operation

1) Running a Recommended Program Assay:

NOTE: It is recommended to allow a five (5) minute rest period between consecutive runs.

1. From the 'Home Screen', select [RECOMMENDED PROGRAMS] [\(Figure V-1\)](#).
2. A list of Recommended Programs has been included in [Appendix 4](#).



Figure V-1: FastPrep-24™ 5G Home Screen

3. Press the appropriate category for the sample type (i.e., plants, bacteria, environmental, human/animal, yeast & fungi, or feces) [\(Figure V-2\)](#).

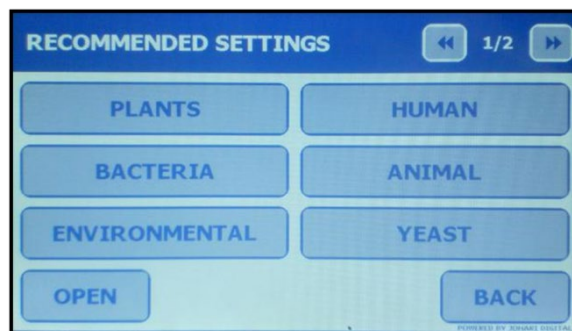




Figure V-2: Recommended Programs, page 1 of 2

4. Scroll through the program listings using the [ / ] buttons located on the top right corner of the 'Touch Screen' [\(Figure V-3\)](#).

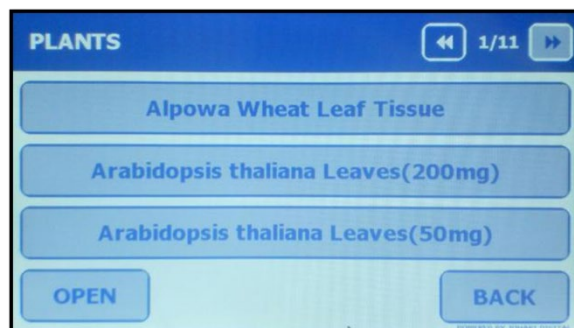


Figure V-3: Plant Category, page 1 of 11

5. Press the desired program button. The FP-24 5G displays the program settings; use the [/] buttons located at the top right corner of the 'Touch Screen' to scroll through each page of parameters. (**Figures [V-4a](#), [V-4b](#) and [V-4c](#)**).

NOTE: Recommended Programs are locked and cannot be edited. Program details are for viewing purposes only.

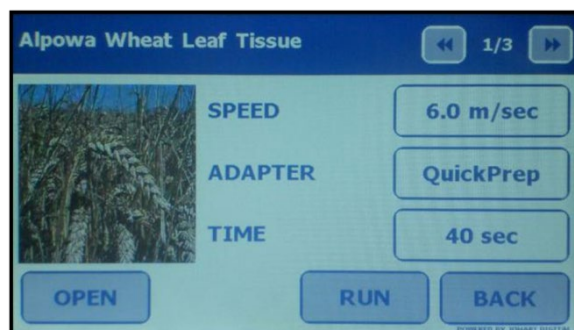


Figure V-4a: Alpowa Wheat Leaf Tissue program, page 1 of 3

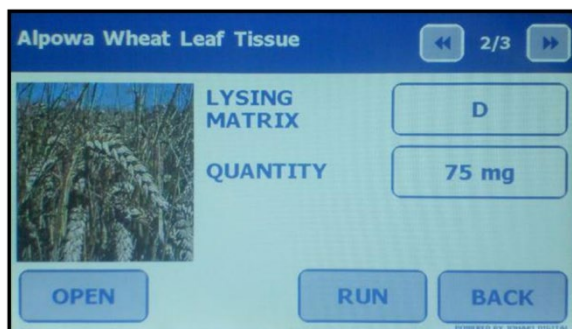


Figure V-4b: Alpow Wheat Leaf Tissue program, page 2 of 3

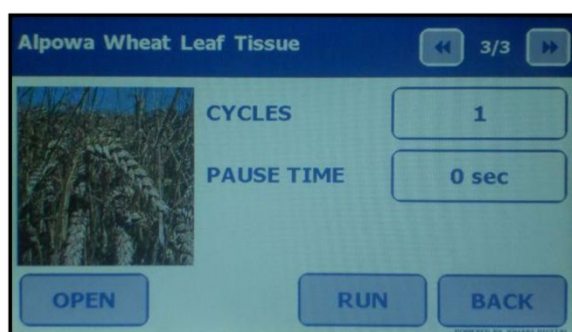


Figure V-4c: Alpow Wheat Leaf Tissue program, page 3 of 3

6. Press [RUN] to initiate the program. The FP-24 5G prompts for confirmation of the program run ([Figure V-5](#)).

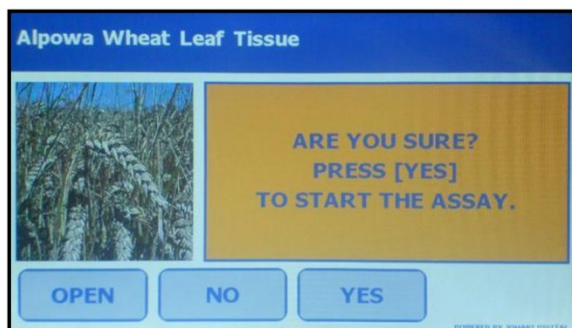


Figure V-5: Confirmation Screen for Recommended Program Assay Run

7. Press [YES] to initiate the assay run. The FP-24 5G starts the assay while displays a countdown of the assay time remaining ([Figure V-6](#)).

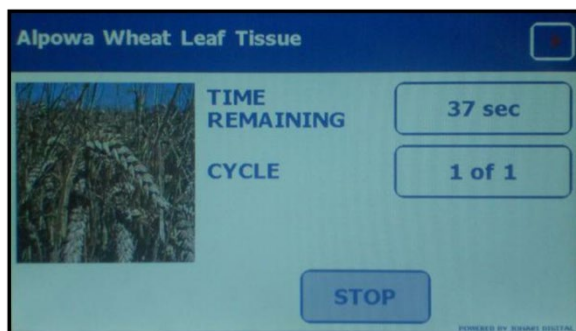


Figure V-6: FastPrep-24™ 5G ‘Touch Screen’ Countdown Display During Assay Run

NOTE: Any assay may be aborted at any time during the assay run by pressing [STOP].

NOTE: If the Program contains multiple Cycles and Pauses, a separate countdown display will show during Pause time ([Figure V-7](#)).

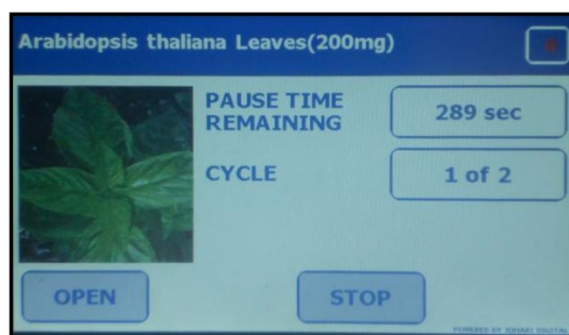


Figure V-7: FastPrep-24™ 5G ‘Touch Screen’ Countdown Display During Cycle Pause

8. Upon completion of the assay, the FP-24 5G screen will display the message “ASSAY COMPLETED SUCCESSFULLY” ([Figure V-8](#)).

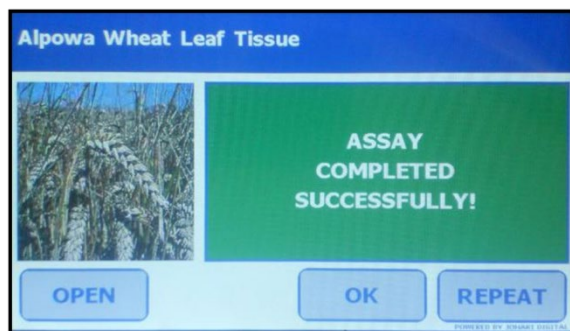


Figure V-8: Confirmation Screen for Completion of Recommended Program Assay Run

9. After an assay has been completed.
 - a. Press [OPEN] to release the dome lock mechanism and retrieve your samples.
 - b. Press [REPEAT] to run the same assay again.
 - c. Press [OK] to return to selected program main screen.

2) Manually Programming an Assay:

The FastPrep-24™ 5G allows the end-user to program customized assays using variable speed, adapter, time, lysing matrix, quantity, quantity unit, cycles, and pause time parameters. To program a custom assay using the FastPrep-24™ 5G:

1. Press [PROGRAM MANUALLY] ([Figure V-1](#)) at the 'Home Screen'. The FP-24 5G Touch Screen will display the first of three total pages of programmable variables ([Figures V-9a](#), [V-9b](#) and [V-9c](#)).

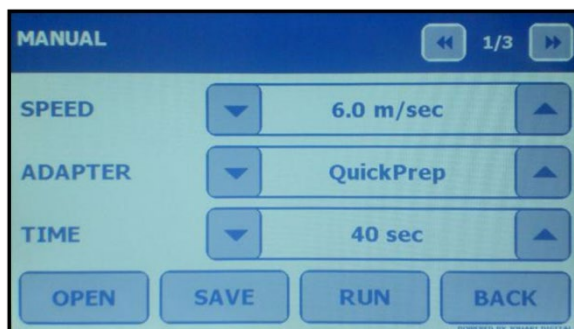


Figure V-9a: Manual Assay Parameters, page 1 of 3

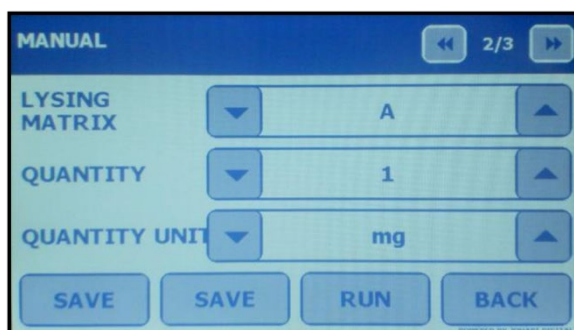


Figure V-9b: Manual Assay Parameters, page 2 of 3

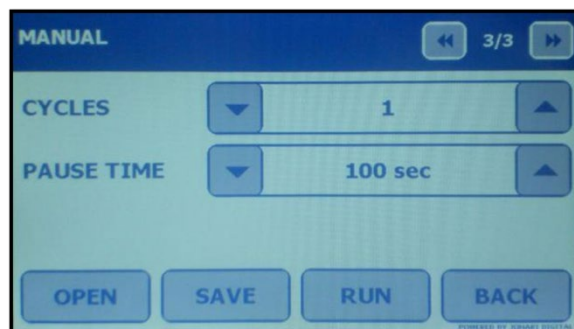






Figure V-9c: Manual Assay Parameters, page 3 of 3

2. Press the up/down arrows [ / ] to program each parameter, using the scroll buttons in the top right corner of the screen to move from page to page. A complete list of parameters, parameter ranges, and default settings is found in the FastPrep-24™ 5G [programmable settings](#) in [Section 1](#).

NOTE: The up/down arrows [ / ] are equipped with a convenient Rapid Scroll Feature. Holding the arrows down will rapidly increase or decrease the desired setting. Pressing and immediately releasing the arrows will increase or decrease by one default setting per action.

NOTE: The Manual Program mode contains three (3) screens of programmable variables. Any assay parameters not programmed will run at default levels.

NOTE: Speed settings higher than 6.5 m/s are available only for the QuickPrep-3 adapters. All other optional adapters are limited to a maximum speed of 6.5 m/s.

IMPORTANT



While the FastPrep-24™ 5G is capable of run times up to 120 seconds, ***it is highly recommended that any run time over 60 seconds be monitored very closely!*** Some Lysing Matrix combinations can generate significant heat after 60 seconds at high speeds and tube failure is possible.

3. Once all parameters are programmed, the end user can choose to save the assay for easy recall or to run the assay.

F. Running an Assay

1. To run the program without saving, press [RUN] to initiate the program. The FP-24 5G prompts “ARE YOU SURE? PRESS [YES] TO START THE ASSAY” for confirmation of the program run ([Figure V-10](#)).

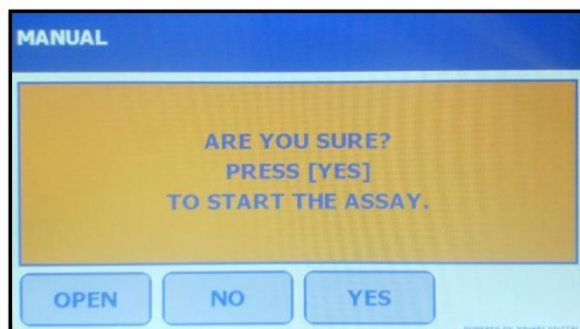


Figure V-10: Confirmation of Program Run

2. Press [YES] to initiate the assay run. The FP-24 5G will start the assay while displaying a countdown of the assay time remaining ([Figure V-11](#)).

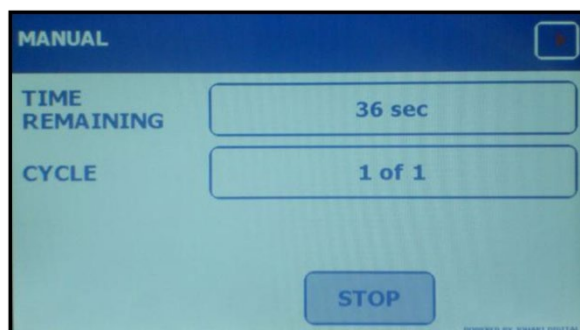


Figure V-11: FastPrep-24™ 5G Countdown Display during Manual Assay Run

3. Upon completion of the assay, the FP-24 5G screen displays “ASSAY COMPLETED SUCCESSFULLY!” ([Figure V-12](#)).

NOTE: All parameters entered in the [Program Manually] mode will be saved until the program is manually edited or the system is turned off. Once power is resumed, settings will revert to factory defaults.

4. After an assay has been completed ([Figure V-12](#)):
 - a. Press [OPEN] to release the dome lock mechanism and retrieve your samples.
 - b. Press [REPEAT] to run the same assay again.

- c. Press [SAVE] to save program.
- d. Press [OK] to return to selected program main screen.

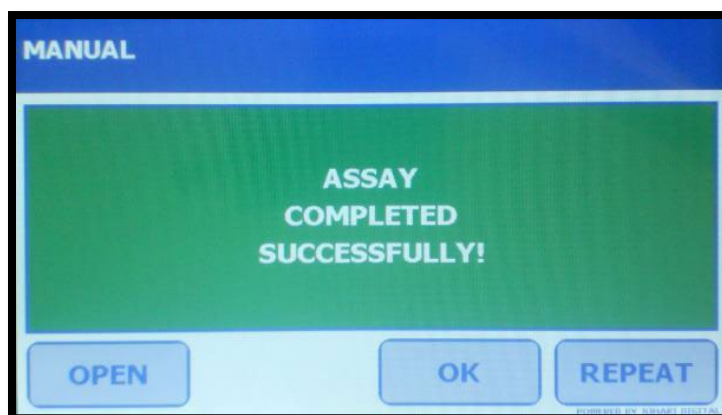


Figure V-12: Manual Assay Complete

G. Saving an Assay

In the Program Manually mode, an assay may be saved immediately after programming ([Figure V-9](#)) or upon completion of the assay ([Figure V-12](#)). To save an assay:

1. Press [SAVE] at the bottom of screen. The FP-24 5G will automatically display a keypad ([Figure V-13](#)).

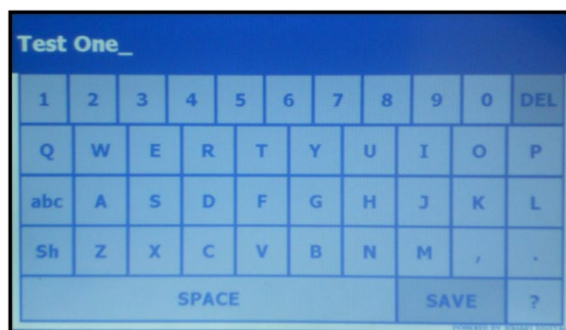


Figure V-13: FastPrep-24™ 5G Keypad

2. Using the keypad, enter the program name. Alpha/numeric and punctuation is limited to 25 characters and spaces.

NOTE: Each saved assay must be given a unique name.

NOTE: The maximum number of saved programs allowable is twelve (12).

3. Press [ENTER] to save named program to memory. The FP-24 5G prompts “ARE YOU SURE YOU WANT TO SAVE {ASSAY NAME}?”. Press [YES] to save.
4. The FP-24 5G displays “{ASSAY NAME} SAVED SUCCESSFULLY!”.
5. Press [OK] to return to the Program Manually screen.
6. To access saved programs again, press [SAVED PROGRAMS] from the FP-24 5G main menu.

NOTE: Each program must be saved under a unique name. Repeating a name will result in a program error “THE ASSAY WITH THE SAME NAME EXISTS. PLEASE CHANGE THE NAME OF ASSAY!”. The name of the last saved program will remain on result line of the keypad entry screen until unit is powered off.



NOTE: A program cannot be edited once saved.

3) Using Saved Programs:

To access the saved programs, press [SAVED PROGRAMS] from the ‘Home Screen’.
[\(Figure V-1\)](#)

A. Retrieving a Saved Program

To retrieve a previously saved program:

1. Press [SAVED PROGRAMS] from the FastPrep-24™ 5G ‘Home Screen’.
2. Scroll through the pages using the   buttons until the program location is found [\(Figure V-14\)](#).

3. Press {ASSAY NAME} to open the program. View the program settings using the scroll [<< / >>] buttons located at the top right corner of the 'Touch Screen'.
4. Press [RUN] to start the program. The FP-24 5G prompts, "ARE YOU SURE? PRESS [YES] TO START THE ASSAY".
5. Press [YES] to start the assay run.
6. When the run is completed, the FP-24 5G displays "ASSAY COMPLETED SUCCESSFULLY!".

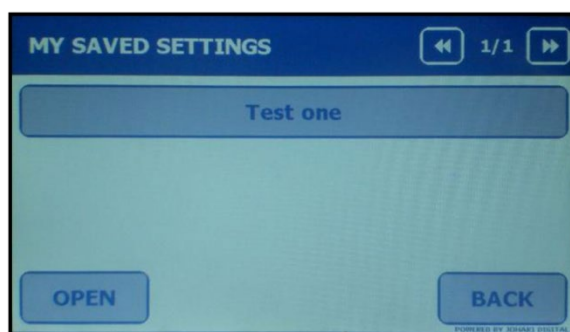


Figure V-14: FastPrep-24™ 5G [SAVED PROGRAMS] Menu Listing

7. After an assay has been completed:
 - a. Press [OPEN] to release the dome lock mechanism and retrieve your samples.
 - b. Press [REPEAT] to run the same assay again.
 - c. Press [OK] to return to selected program main screen.

NOTE: A program cannot be edited once saved.

B. Deleting a User-Programmed Assay:

1. Press [SAVED PROGRAMS] from the 'Home Screen' ([Figure V-1](#)). Scroll through the pages using the scroll [<< / >>] buttons located at the top right corner of the 'Touch Screen'.
2. Press the {ASSAY NAME} to be deleted.

3. Press [DELETE] to delete program. The FP-24 5G displays “ARE YOU SURE YOU WANT TO DELETE {ASSAY NAME}” to confirm deletion. Press [YES] to delete. [\(Figure V-15\)](#)

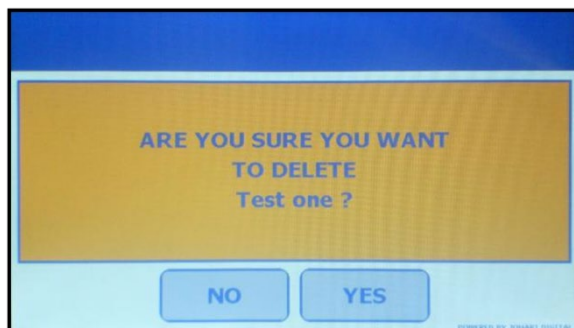


Figure V-15: Confirm Delete Program (Assay Name “Test one”)

4. Once the program is deleted, the FP-24 5G displays “{ASSAY NAME} DELETED SUCCESSFULLY!”. [\(Figure V-16\)](#)

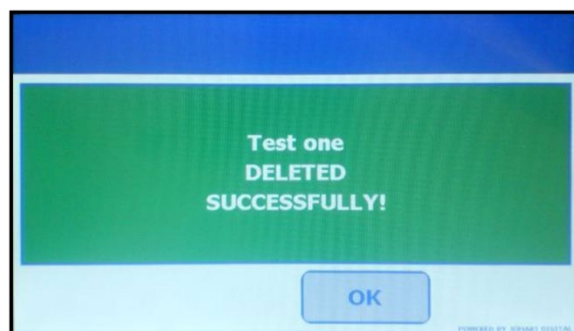


Figure V-16: Program Deleted Successfully (Assay Name “Test one”)

5. Press [OK] to return to previous screen.

4) **Stopping Cycle in Progress:**

A cycle in progress on the FP-24 5G can be stopped as needed. To stop a cycle in progress:

1. Press the [STOP] button located at the bottom of the screen [\(Figure V-11\)](#). The FP-24 5G prompts “ARE YOU SURE YOU WANT TO ABORT ASSAY” [\(Figure V-18\)](#). Press the [YES] button to abort the assay, or press [NO] to return to the previous screen.

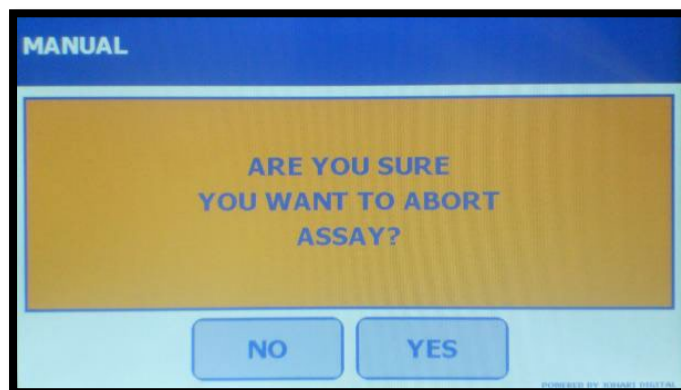


Figure V-18: Assay Abort Screen

2. If an assay is aborted, the FP-24 5G displays a confirmation screen, prompting “SESSION IS ABORTED!” [\(Figure V-19\)](#). Press [OK] to return to the main menu or [OPEN] to open the Dome.

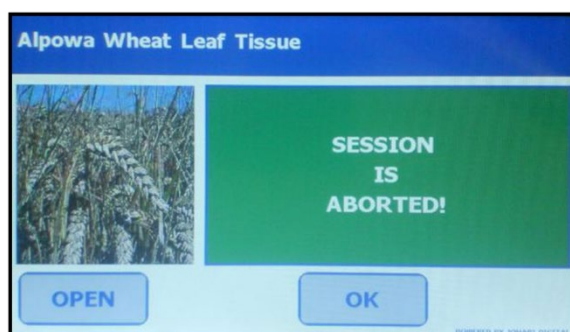


Figure V-19: Confirmation Screen for Aborted Assay

B. FastPrep-24™ 5G Administrative Functions

The FastPrep-24™ 5G allows end users to activate or deactivate system features, export data records, install software upgrade and access other onboard information. To access Administrative Functions [\(Figure V-20\)](#), press [ADMIN TOOLS] on the ‘Home Screen’

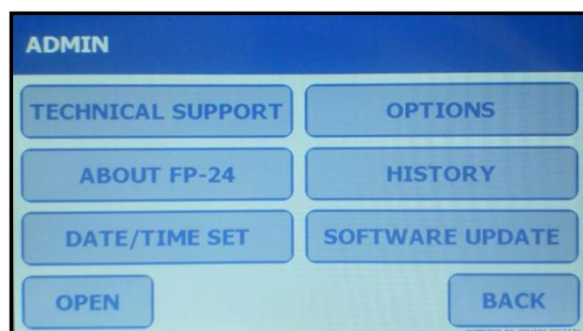


Figure V-20: [ADMIN TOOLS] Main Menu

1) **About FastPrep-24™ 5G: System Versions and End-User Agreement**

The 'About FP24' screen provides information on the FastPrep-24™ 5G, such as current software and firmware versions, as well as the end user agreement. To access information:

1. Press the [ABOUT FP24] button. The FP-24 5G will display the current application (software) and driver (firmware) versions [\(Figure V-21\)](#).

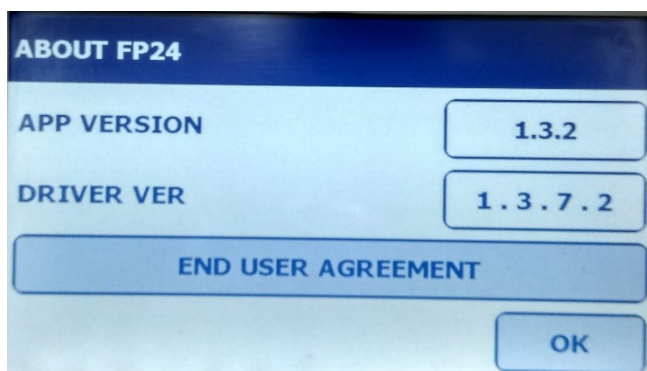


Figure V-21: About FastPrep-24™ 5G

2. Press [END USER AGREEMENT] to view the MP Biomedicals End User Agreement [\(Figure V-22\)](#). Press the [UP]/[DOWN] buttons to scroll through the text.
3. Read and accept the terms of the End User Agreement by pressing [OK]. This will exit the 'End User Agreement' Screen and return to the About FP-24 screen.

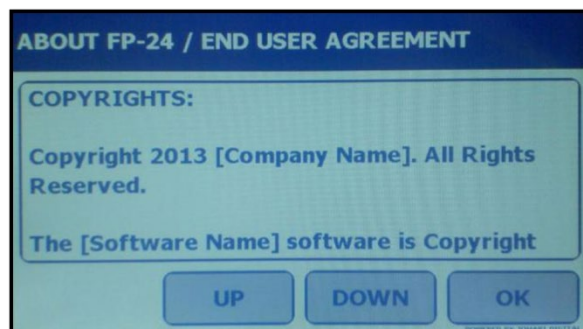


Figure V-22: End User Agreement

4. Press [OK] to return to the 'Admin Tools' menu.



NOTE:

After accepting the terms, please register your New FastPrep-24™ 5G at

<https://www.mpbio.com/us/registermyfastprep>

2) Date /Time Set

The 'Date / Time Set' menu allows the end-user to set the date and time on the instrument. The date is formatted using month, day and year, and the time is displayed using a 24 hour clock.

1. Press the [DATE/TIME SET] button on the 'Admin Tools' main menu.
2. Press the up/down arrows [ / ] to program each parameter [\(Figure V-23\)](#).

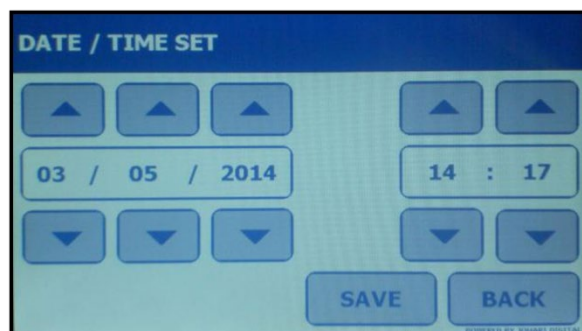


Figure V-23: 'DATE/TIME SET' Screen

3. Press [SAVE] to save changes. The FP-24 5G prompts "ARE YOU SURE YOU WANT TO SAVE NEW PARAMETERS?". Press [YES] to save. [\(Figure V-24\)](#).

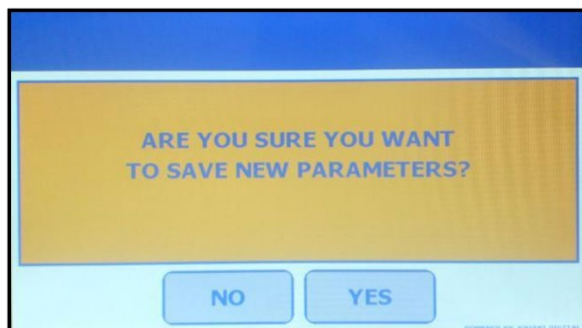


Figure V-24: Confirm Parameters Screen

4. The FP-24 5G will prompt “NEW PARAMETERS SAVED SUCCESSFULLY” ([Figure V-25](#)). Press [OK] to return to ‘Admin Tools’ menu.

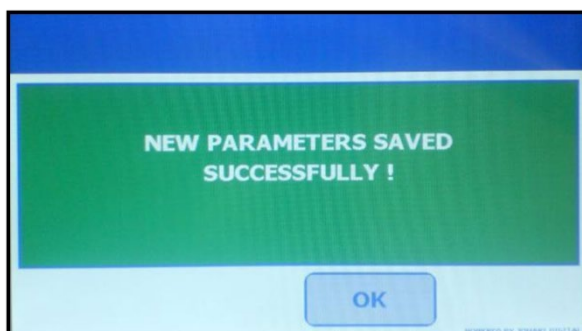




Figure V-25: New Parameters Saved Successfully Screen

3) Optional Settings

Additional FP-24 5G features, including the strobe light, keyboard clicks and voice interface, may be accessed through the ‘Options’ menu. To access the additional FP-24 features:

1. Press [OPTIONS] on the [ADMIN TOOLS] main menu.
2. Scroll through the pages using the [ / >] buttons located at the top right corner of the ‘Touch Screen’ ([Figure V-26a](#) and [V-26b](#)).

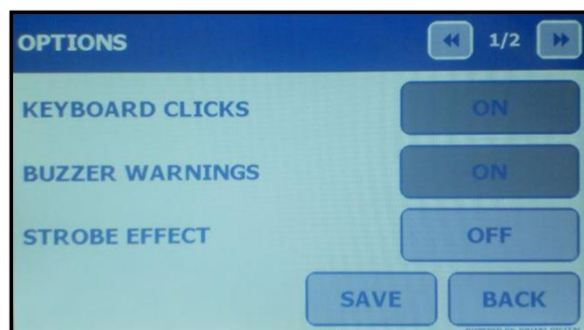


Figure V-26a: Options Screen (Page 1)

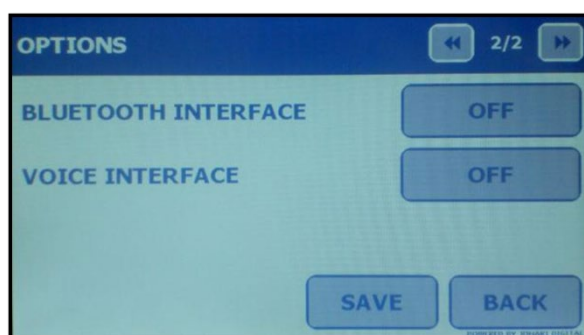


Figure V-26b: Options Screen (Page 2)

A. Keyboard Clicks

The 'Keyboard Clicks' feature controls the level of sound made when a button is pressed on the 'Touch Screen'. The Keyboard Clicks feature is controlled using a toggle button; press [ON] or [OFF] to turn on the sound or to silence respectively.

B. Buzzer Warnings

The 'Buzzer Warnings' feature controls the level of sound made for various minor confirmatory indication or out of compliance parameters. These **warnings** always accompany a **Yellow Warning Screen** and the user has the option to continue/over-ride. The 'Buzzer Warnings' feature is controlled using an [ON/OFF] toggle button: [ON] activates while [OFF] silences this feature.

Examples of these **warnings** include:

Program parameter is not acceptable.

Assay aborted in session

New administrative parameters saved	Assay deleted
Assay initiated	Data storage near capacity
Assay time complete	FastPrep-24™ 5G instrument warm
FastPrep-24™ 5G overloaded (current)	



More serious Audio System **Alarms** are accompanied by a **Red Alarm Screen**; The user cannot over-ride these, and these cannot be deactivated. Example: System Overheating.

C. Strobe Effect

The 'Strobe Effect' feature controls the two strobe lights located within the sample chamber. The 'Strobe Effect' is useful for optimizing assay parameters because it allows for visual inspection of sample lysis progression in real time. The 'Strobe Effect' feature is controlled using a toggle button. Press [ON] to enable strobe lighting and [OFF] to disable this feature.

NOTE: Bluetooth Interface and Voice Interface Options are currently unavailable at this time and are therefore disabled. Once available, these icons will be activated.

4) Software Update

The FastPrep-24™ 5G features custom software updates. Registered end-users receive software updates via email or portable drives, and the FastPrep-24™ 5G software can then be updated using a USB port directly located directly on the instrument along with a standard laptop installed with the FastPrep-24™ 5G Host Application.

Materials Required

- FastPrep-24™ 5G instrument
- Laptop installed with FastPrep-24™ 5G Host Application
- USB A/Mini B Cable (provided)
- Pin, such as push pin, straightened paper clip, or other thin, rigid wire

NOTE: The FastPrep-24™ 5G Host Application is currently compatible with MS Windows 10 or MS Windows XP operating systems. Contact MP Bio Technical Service for other operating systems.

a) Installing the FastPrep-24 5G Host Application

Registered end-users will receive application files for the FastPrep-24 5G Host Application as zipped files via email or portable drive.

1. Download the attached zip file onto the computer.
2. Unzip the setup.exe and .dfu files on the desktop of the computer.
3. Install the program on the computer by double click on the setup.exe file and follow the instructions on the new window.
4. Shortcut icon for program will be created on the desktop, called “FastPrep-24 5G” (Figure V-27).



Figure V-27 : FastPrep-24 5G Program icon

5. Program is now installed

b) Install FastPrep-24 5G driver

1. Double click the FastPrep-24 5G Program icon (Figure V-27).
2. FastPrep-24 5G program window will open.
3. Next, on the desktop, right click on the FastPrep-24 5G Program icon (Figure V-27) and select “Properties” from the dropdown window (Figure V-28).

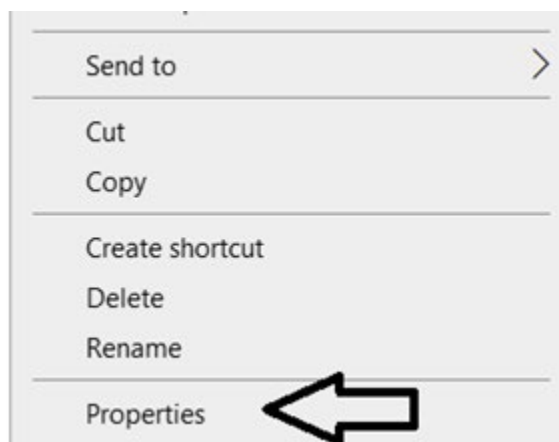


Figure V-28: Dropdown window to select Properties

4. On the new window (Figure V-29), under the tab “Shortcuts,” select “Open File Location.”

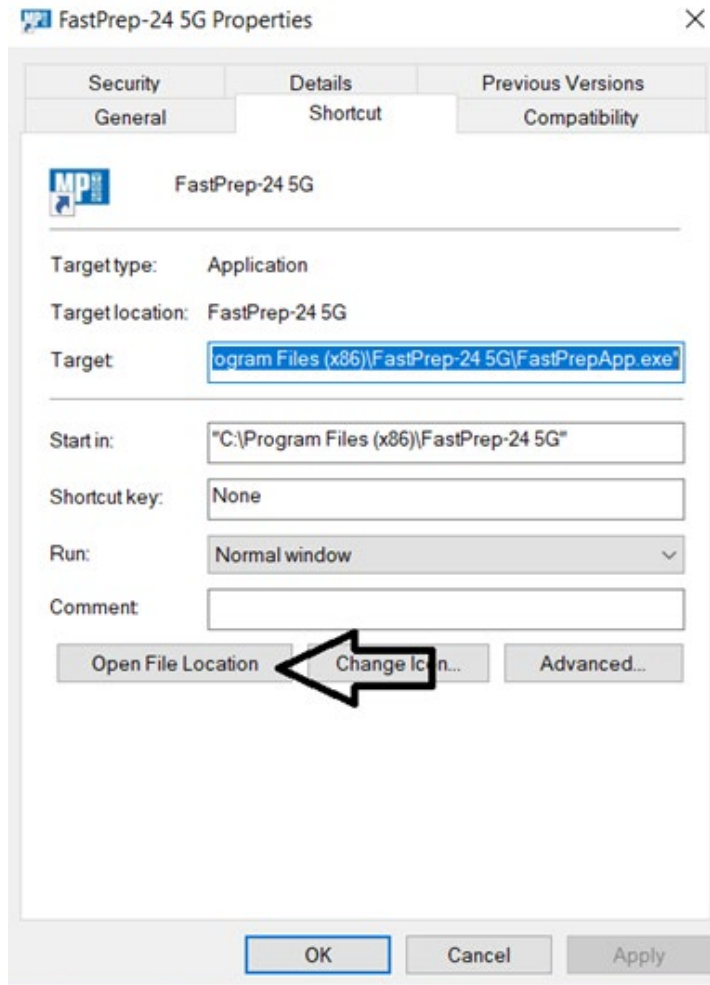


Figure V-29: FastPrep-24 5G Properties Window

5. On the new dialog box window (Figure V-30a), select the Driver folder and find the application files.

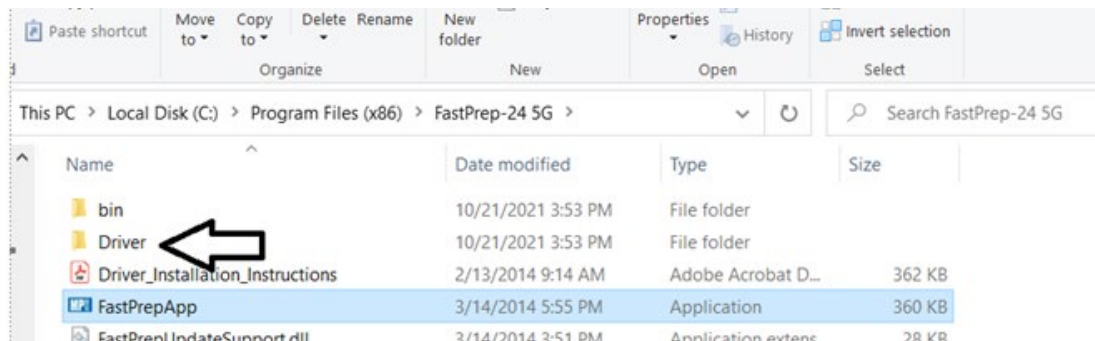


Figure V-30a: Dialog Box Window

- Right click on the STtube file and from the dropdown window select Install (Figure V-30b).

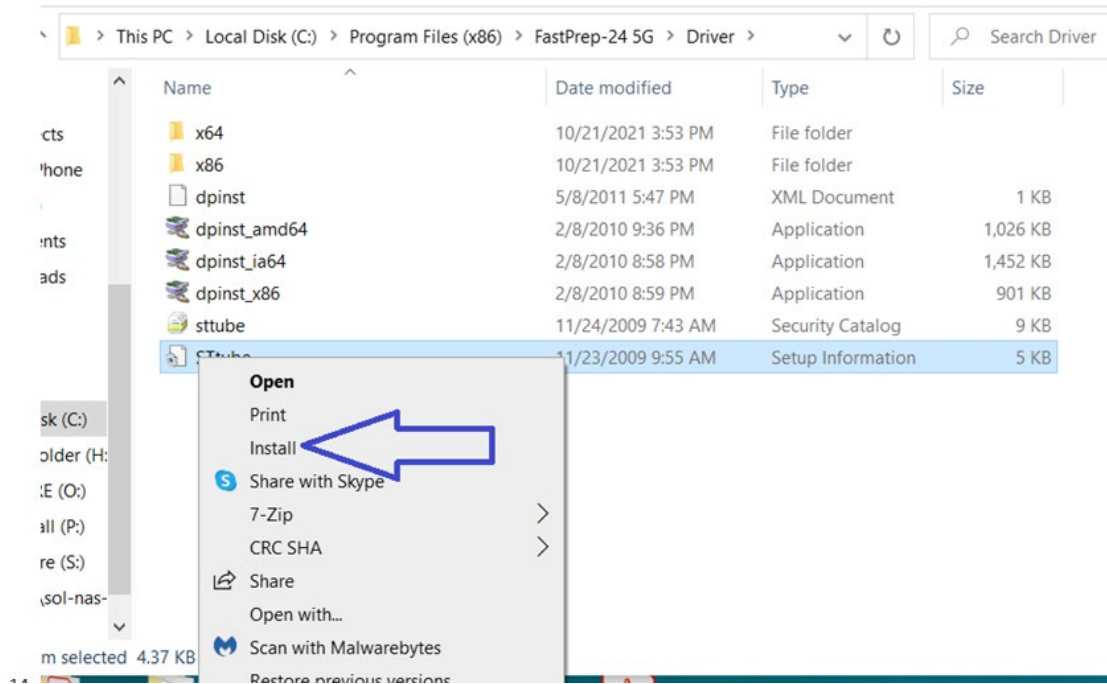


Figure V-30b: Dialog Box Window for the installing application file

- Window will display the message “The operation completed successfully.” Driver is now installed.

c) Open the FastPrep-24 5G in Safe Mode

1. FastPrep-24 5G unit should be turned Off when starting.
2. As the FastPrep-24 5G is not connected, the Host Application will not detect a device and no buttons [Update Firmware] and [Read Report from Device] will be activated (Figure V-31).



Figure V-31: FastPrep App Home Screen (device not detected).

3. Connect the USB cable first to the laptop using the large Type A plug, then to the FP-24 5G using the smaller Type B plug.
4. Depress the “Safe Mode Boot” button adjacent to the USB Port (Figure V-32). Insert safety pin to depress this button while turning the unit to ON. The FastPrep-24 5G will start up in Safe Mode.

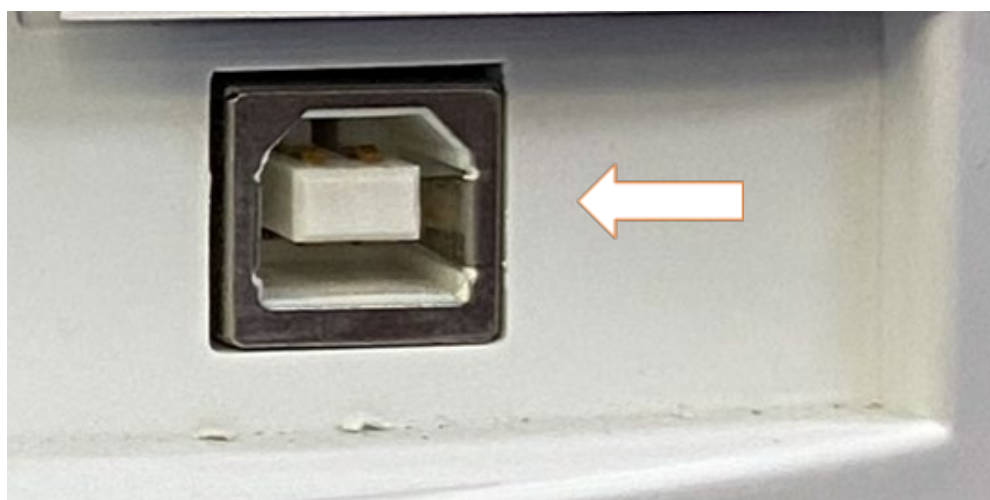


Figure V-32: “Safe Mode Boot” Button found on FastPrep-24 5G Rear Panel

5. Depress the 'Safe Mode Boot Button' until the FastPrep-24 5G Host Application detects the attached device. This can take several seconds. The Update Firmware will be activated, and the FastPrep-24 5G LCD screen will be dark. Once device detection occurs, release the 'Safe Mode Boot Button' by removing the pin.

d) Install the Firmware Update

1. Press the activated button [Update Firmware] on the FastPrep-24 5G Touch Screen.
2. On the new Dialog Box window, select the binary file (.dfu) found on the desktop and click [OPEN] to start installation.
3. It will take about 2 minutes to go through the steps of Erasing, Upgrading and Verifying (Figures V32a, V-32b and V32c).



Figure V-32a: FastPrep App Progress Bar (Erasing)

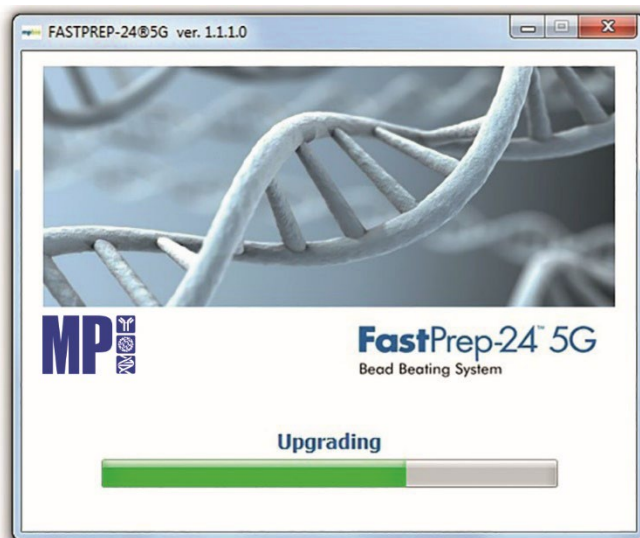


Figure V-32b: FastPrep App Progress Bar (Upgrading)



Figure V-32c: FastPrep App Progress Bar (Verifying)

4. The FP-24 5G Host Application displays a Dialog Box (Figure V-33) upon successful completion of the file upload. Click [OK] to exit.



Figure V-33: FastPrep App Update Successfully Completed

5. Reboot the FastPrep-24 5G unit by powering off for a minimum of 30 seconds before powering back on.
6. Power on the FP-24 5G in normal (standard) mode.
7. FastPrep-24 5G will launch upgrade software progress bar on the LCD screen.
8. Upon completion, the screen will display “Drive MCU Firmware Updated Successfully”. Press [OK].
9. Reboot the FastPrep-24 5G unit by powering off for a minimum of 30 seconds before powering back on.
10. Software Firmware is now updated.
11. Verify a successful upgrade installation using the [ABOUT FP24] button found under [ADMIN TOOLS]. The ‘App Version’ and ‘Driver Version’ will correspond to the new upgrade([Figure V-34](#)).
12. Additional verification and validation activities may be required. MP Biomedicals strongly recommends testing using the following functions before running samples: running an assay, saving an assay, opening the dome using the Touch Screen, and changing an optional setting.

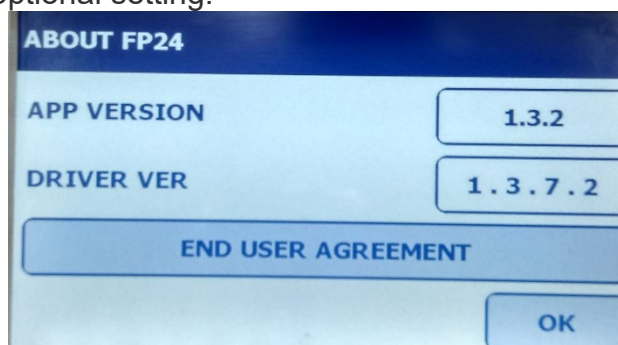


Figure V-34: About FastPrep-24 5G

e) Calibrating the FastPrep-24™ 5G ‘Touch Screen

The ‘Touch Screen’ may require calibration upon completion of software update. In cases requiring calibration, the ‘Touch Screen’ prompts, “Touch Panel Calibration- Step No. 1 of 4.” Follow the instructions on each screen, pressing the locations as indicated to calibrate touch screen” ([Figure V-35](#)).

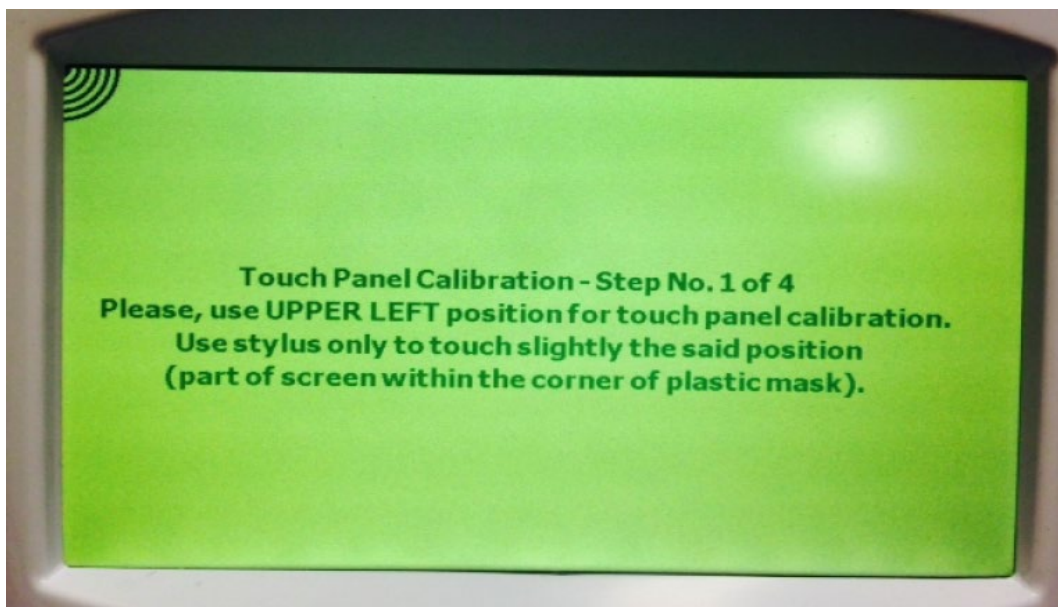


Figure V-35: FastPrep-24™ 5G ‘Touch Screen’ Calibration Prompt

To calibrate the ‘Touch Screen’, use a stylus or pencil to firmly press the locations as indicated. The positions will appear in sequence in each quadrant, beginning with the upper left corner ([Figure V-35](#)), followed by the lower left corner, the lower right corner and finally the upper right corner.

NOTE: Use firm but not hard force when depressing the stylus or pencil during ‘Touch Screen’ calibration. Use of too much force can damage the ‘Touch Screen’.

5) Exporting History File

The FastPrep-24™ 5G stores assay information that is available for export as needed to a laptop or PC. Historical data is exported in a CSV format that contains information such as assay date and time, assay name and category, assay parameters (e.g., speed, time, cycles, etc.) and whether the assay completed or aborted. While the FastPrep-24™ 5G can store information on up to 2048 assay, it is recommended that the system history be exported and/or purged routinely. To export an Assay History File:

1. Launch the FastPrep-24™ 5G Host Application by double clicking the FastPrepApp.exe program file or the program icon on your desktop.

NOTE: If the FastPrep App has not been installed on the laptop/PC, please refer to Section 4 above, UPDATING SOFTWARE VERSIONS, Steps 1-3 for application installation.

2. Power on the FP-24 5G at the 'Rocker Switch' [\(Figure I-2\)](#).
3. Connect the USB cable first to the laptop using the large Type A plug, then to the FP-24 5G using the smaller Type B plug [\(Figure V-29\)](#).
4. After a few seconds, the FastPrep App detects the device and activates the [READ REPORT FROM DEVICE] button [\(Figure V-36\)](#).



Figure V-36: FastPrep App Showing Device Detected

5. Click [READ REPORT FROM DEVICE] to initiate file download. A Dialog Box will open allowing you to select a location to save the CSV file.
6. Rename the file as appropriate and select the desired save location. Click [SAVE] to save.
7. The FP-24 5G Host Application then prompts for the purging of stored data from the FP-24 5G by displaying the message "Do you want to delete all entries in history?" [\(Figure V-37\)](#). Press [YES] to clear the data history cache or [NO] to continue storing data.

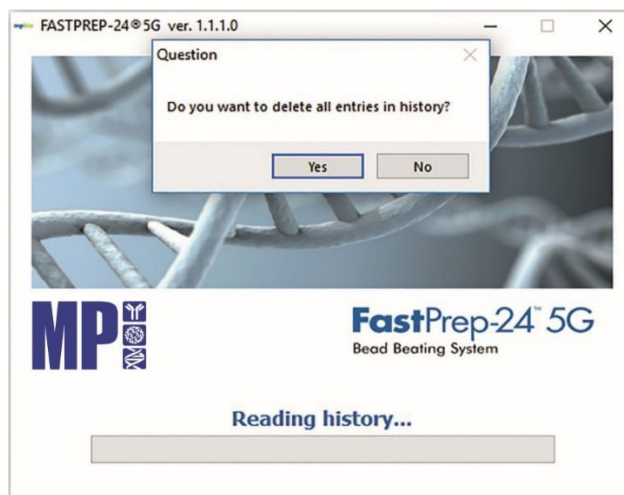


Figure V-37: Purge History using FastPrep-24™ 5G Host Application

8. Alternatively, to clear the data storage from the FP-24 5G without previously saving, press [ADMIN TOOLS] from the main menu, followed by [HISTORY]. On the History Screen, press [ERASE ALL RECORDS] to purge the data file ([Figure V-38](#)).

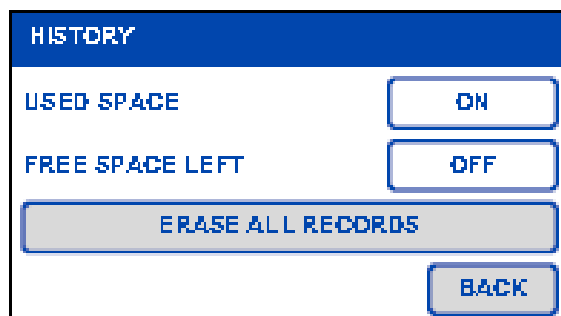


Figure V-38: Purge File from History Screen

9. The FP-24 5G prompts for “ARE YOU SURE YOU WANT TO DELETE HISTORY LOG?”. Press [YES] to erase ([Figure V-39](#)).

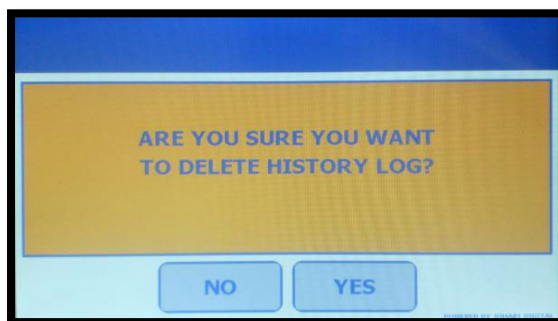


Figure V-39: Purge Confirmation Screen

10. The FP-24 5G prompts for "HISTORY LOG DELETED SUCCESSFULLY?". Press [OK] to return to the previous screen ([Figure V-40](#)).

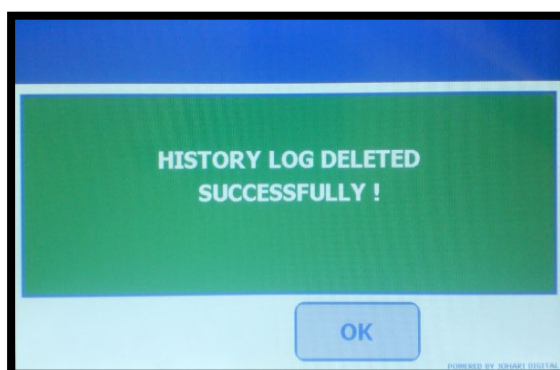


Figure V-40: Purge Successful Screen

6) Technical Support

The FastPrep-24™ 5G contains onboard contact information for all of MP Biomedicals' technical services on a global basis; regions are listed by country. These technical services support all aspects of your FastPrep instrument, including technical support and supply of FastPrep-24™ 5G consumables such as Lysing Matrix Tubs and Purification kits. To access 'Technical Support':



1. Press [TECHNICAL SERVICE] from the 'Admin Tools' menu ([Figure V-21](#)).
2. Scroll through the different regions listed using the scroll buttons [ / ] located at the top of the screen ([Figure V-41](#)).



Figure V-41: Global Technical Services Listing by Country (page 1 of 4 shown)

3. Press [COUNTRY NAME] to display the desired MP Biomedicals' Technical Services contact information, including email address and telephone number ([Figure V-42](#)).

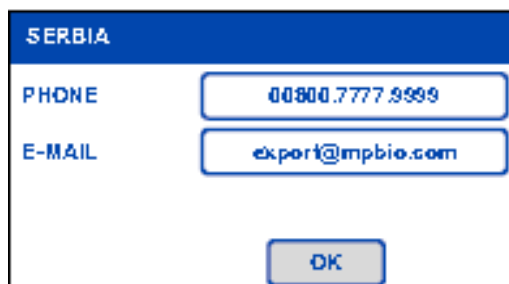


Figure V-42: Country Specific Service Listing

Resetting the FastPrep-24™ 5G:

The FastPrep-24™ 5G is designed to retain all saved programs upon loss of power to the instrument (e.g., emergency stop deployed, unplugging the instrument, etc.). The instrument will revert to default settings for the manual programming mode.

SECTION V: WARRANTY & LIABILITY

The FastPrep-24™ 5G is warranted against defects in material and workmanship for one (1) year after the date of delivery to the original purchaser. This warranty is limited to defects in materials and workmanship and does not cover incidental or consequential damages.

MP Biomedicals, will repair, free of charge, any apparatus covered by this warranty. This warranty includes one-year parts and labor in MP Biomedicals' facilities or by approved distributors. Warranty work is subject to the inspection of unit. No instruments, equipment, or accessories will be accepted without a Return Material Authorization (RMA) number issued by MP Biomedicals. Costs of shipping the unit are not covered in this warranty. The warranty obliges you to follow all precautions as listed in this manual; failure to do so will void warranty.

An instrument that may contain hazardous and / or infectious materials must be packed and labeled according to the U.S. Department of Transportation (DOT) and / or European Community (EC) regulations applying to the transportation of hazardous and / or infectious materials. All shipping documents must meet DOT and / or EC regulations. All returned units must be fully decontaminated of any chemical, biological or infectious agents.

Use of the FastPrep-24™ 5G in a manner other than that specified in this user manual may jeopardize personal safety. MP Biomedicals shall not be liable for damages due to the improper handling, abuse, or unauthorized repair of this instrument. MP Biomedicals, LLC assumes no liability, express or implied, for use of this instrument.

MP Biomedicals supplies kits, reagents and adapters for use with the FastPrep-24™ 5G. Use of kits, reagents and Adaptors not supplied by MP Biomedicals with the FastPrep-24™ 5G is not covered under this warranty.

SECTION VI: MAINTENANCE

The FastPrep-24™ 5G is a robust instrument. Depending on the intensity and frequency of use certain wear parts and moving parts must be checked and are likely to require a long-term replacement to keep the FastPrep-24™ 5G in good working condition.

These are parts that undergo movement or friction.

Contact MP Biomedicals in the event of an instrument question or to subscribe for a yearly maintenance contract.

- An annual revision of your instrument
- Assistance within 48 hours
- An instrument always available
- A service tailored to your own needs

A. Maintenance

The FastPrep-24™ 5G should be cleaned following use of the instrument, or if a spill occurs during use. If a sample tube leaks during a cycle, the inside of the 'Dome' and surrounding areas may be contaminated. To clean the FastPrep-24™ 5G:

1. Turn the FP-24 5G off and disconnect the power cord from the instrument.
2. Wipe the exterior of the FP-24 5G with a paper towel moistened with 7X™ Cleaning Solution.
3. Dry the exterior of the FP-24 5G with a dry paper towel.
4. Open the dome and remove the Adaptor.
5. Remove the Bowl.
6. Clean the Adaptor and the bowl using 7X™ Cleaning Solution, rinse with distilled water, and dry using paper towels.
7. Wipe the interior of the FP-24 5G with a paper towel moistened with 7X™ Cleaning Solution.
8. Wipe the interior of the FP-24 5G with a damp paper towel to remove any residual 7X™ Cleaning Solution.
9. Dry the interior of the FP-24 5G with a dry paper towel.
10. Replace the Adaptor and the bowl than close the dome
11. Regularly clean the device including area under the bowl.

DAMAGE TO INSTRUMENT

Use of a cleaning agent other than 7X™ Cleaning Solution may damage the FastPrep-24™ 5G. Bleach, solvents, or acidic, alkali or abrasive reagents may damage the gel coat finish, cause corrosion of metal parts, or crazing of the transparent lid.

7X™ Cleaning Solution

MP Biomedicals recommends that FastPrep™-24 5G cleaning be performed using the MP Biomedicals 7X™ Cleaning Solution. 7X™ Cleaning Solution has been a safe and effective cleaning solution for many laboratory applications.

7X™ Cleaning Solution is composed of a proprietary mixture of a strong, phosphate-free sequestering reagent, a completely soluble and non-toxic solvent, and powerful, environmentally friendly wetting / foaming / emulsifying agents. This combination, along with powerful anionic surfactants, makes 7X™ Cleaning Solution an ideal solution for applications that require low cell toxicity, low debris, high efficiency cleaning, and no corrosion. As such, 7X™ Cleaning Solution is ideal for thorough cleaning of sensitive equipment such as, but not limited to the FastPrep-5G, bioreactors, culture apparatus, and general laboratory equipment including test tubes, microscopy slides, pathology equipment, burets, pipettes, etc.

Regular 7X™

With extremely powerful and effective surfactants, this powerful solution was designed to clean laboratory plastics and glassware by simply soaking overnight. Alternately, equipment may be boiled.

Item	Catalog No.
1 Gallon Plastic Bottle	097667093
4 x 1 Gallon Case	097667094

ES-7X™

“Environmentally Safe” phosphate-free ES works just like the original 7X™ and is designed to work in soaking or non –agitated applications. ES-7X™ is highly concentrated and a little goes a long way.

Item	Catalog No.
1 Gallon Plastic Bottle	097667193
4 x 1 Gallon Case	097667194

Regular 7X™ Ready-to-Use

A 5% solution of the classic 7X™, pre-diluted for a convenient working solution. Includes a hand pump for easy dispensing.

Item	Catalog No.
1 Gallon Plastic Bottle	097668093

ES-7X™ Ready-to-Use

A 5% solution of the ES-7X™, pre-diluted for a convenient working solution. Includes a hand pump for easy dispensing.

Item	Catalog No.
1 Gallon Plastic Bottle	097668193

B. Fuse Replacement (as needed)

The FastPrep-24™ 5G fuse will require replacement occasionally.

**IMPORTANT**

Disconnect input power cord behind instrument before replacing fuse.

**IMPORTANT**

Use only the specified type of fuse for fuse replacement. Use of non-specified fuses may increase fire danger.

To replace a blown fuse:

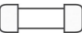
1. Turn the FastPrep-24™ 5G off and unplug the instrument.
2. The 'Fuse Holder' is located directly below the 'Power Inlet Receptacle' and is marked with the  symbol. To remove insert a thin-blade screwdriver or similar into the access hole and gently pry outward, as seen in [Figure VI-1](#).



Figure VI-1: FastPrep™-24 5G Fuse Assembly

3. Remove the faulty fuse and replace it with a new one.
4. Insert the fuse assembly back into the AC receptacle carefully, ensuring that the voltage is correctly configured.
5. Reconnect the power cord and power on the FP-24 5G.

NOTE:


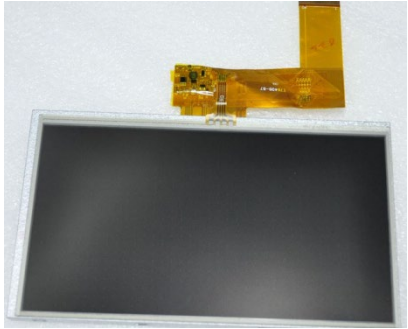



Two (2) spare fuses are provided in the fuse assembly. Two (2) additional fuses are provided with the FastPrep-24™ 5G as part of the accessories. Additionally, fuses may be ordered as needed.

SECTION VII: PARTS INFORMATION





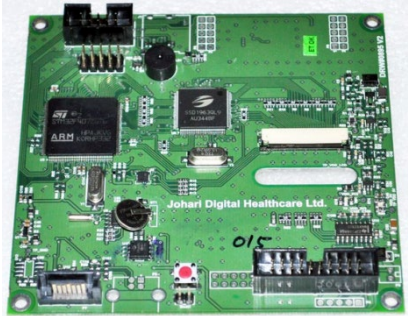



The following spare parts can be ordered directly from MP Biomedicals,






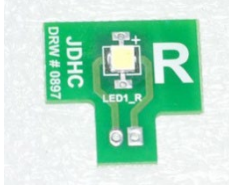



Parts must be replaced by a trained technician.

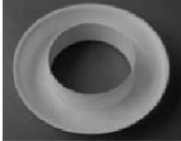Part Description	Catalog Number	Picture
LCD Screen and Touchpad Set 	116005560	
Front Panel Molding	116005565	
Digital (Mother) Board (PCB) 	116005561	

Part Description	Catalog Number	Picture
<p>Driver Board (PCB)</p> 	<p>116005503</p>	
<p>SMPS 48V</p> 	<p>116005502</p>	
<p>Motor with Shaft</p> 	<p>116005504S</p>	

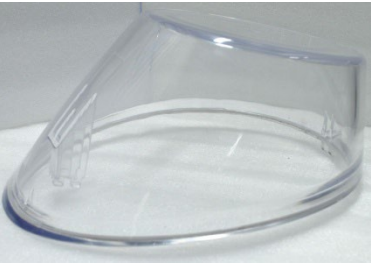Part Description	Catalog Number	Picture
Emergency Stop 	116005524	
Emergency Stop Molding	116005524A	
Strobe Light PCB/Assembled 	116005568: 116005568L Enclosure plastic slight cap Left 116005568R Enclosure plastic slight cap Right	
USB Interface PCB 	116005569	
Bowl – 212 mm	116005505	

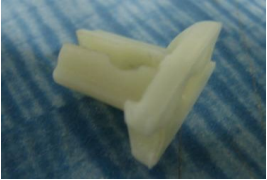

Part Description	Catalog Number	Picture
Electromechanical Latch 	116005563	
Polycarbonate-Dome with Stopper	116005506: Units from 2014 to 2017 116005516 Units 2018 and beyond	
Dome Rubber Gasket	116005507	
Striker Bolt	116005508	
Manual Dome Release Cable Assembly	116005567	
Dome Lock	116005509	

Part Description	Catalog Number	Picture
Dome Rest Assembly Kit	116005589N	 <p>The image shows the components of a dome rest assembly kit. On the left, there is a white plastic bracket, a metal bolt, a white plastic cap, and a black rubber pad. On the right, there is a complete assembly of these parts mounted on a blue base.</p>
Cam Lock	116005588	 <p>The image shows a silver metal cam lock with a handle and a keyhole.</p>
Ratchet Bolt (Length = 25 mm)	116005511	 <p>The image shows a brass ratchet bolt with a threaded shaft and a circular base with two holes.</p>
QuickPrep-3 Adaptor	116005512	 <p>The image shows a white plastic adaptor with a central hole and many small holes around the perimeter.</p>

Part Description	Catalog Number	Picture
<p>3 STEP-Assembly with studded bearings</p> 	<p>116002513</p>	
<p>Spring Assembly Set with Hooks</p>	<p>116002514</p>	
<p>Rubber Shoe</p>	<p>116005515</p>	
<p>Fuse, 10 Amp-UL</p>	<p>116002516</p>	
<p>European AC Cord</p>	<p>1130000062</p>	
<p>3 Pin Flat AC Cord- USA UL</p>	<p>116002519</p>	

Part Description	Catalog Number	Picture
AC Receptacle with Fuse Chamber	116002520	
Rocker Switch-UL	116002521	
Rear Panel Molding	116005570	
Gasket (OD = 150 mm, ID = 45 mm)	116004535	
Inner Ring	116004541	
Outer Ring	116004542	

MAINTENANCE



Part Description	Catalog Number	Picture
Shroud Set	116005564	
USB Cable	116005587	

APPENDIX 1: FASTPREP® INSTRUMENTS

APPENDIX 1: FASTPREP® ADAPTORS

LARGE-SCALE ADAPTORS

Our large sample volume Adaptors are ideally suited for DNA and RNA isolation, enzyme isolation and protein production, natural products isolation, food preparation for quality analysis, biopharma manufacturing, and forensic applications. A wide range of disposable 4.5 ml, 15 mL and 50 mL Lysing Matrix tubes ensure thorough homogenization of any sample type in seconds.



<p>BigPrep™ 2 x 50mL Adaptor Large-scale Adaptor holds 2 x 50 mL Lysing Matrix tubes (compatible with SafTest™ Food Inspection System).</p>	
<p>TeenPrep™ 12 x 15mL Adaptor Medium-size Adaptor holds 12 x 15 mL Lysing Matrix tubes (compatible with SafTest™ Food Inspection System).</p>	
<p>TallPrep™ 24 x 4.5mL Adaptor Large-scale adaptor holds 24 x 4.5ml Lysing Matrix tubes.</p>	

Adaptor Ordering Information			
Cat No	Item	Unit	Price
116002525	BigPrep™ 2 x 50mL Adaptor	1 ea	Contact for details
116002526	TeenPrep™ 12 x 15mL Adaptor	1 ea	Contact for details
116002540	TallPrep™ 24 x 4.5mL Adaptor	1 ea	Contact for details

APPENDIX 1: FASTPREP® INSTRUMENTS

HIGH THROUGHPUT ADAPTORS

Our high-throughput Adaptors are ideally suited for high-throughput applications, up to 24 and 48 samples simultaneously.

<p>HiPrep™ 48 x 2mL Adaptor High-Throughput Adapter holds 48 x 2ml Lysing Matrix tubes.</p>	
<p>QuickPrep-1™ 24 x 2 mL Adaptor</p>	

Adaptor Ordering Information			
Cat No	Item	Unit	Price
116002527	HiPrep™ 48 x 2mL Adaptor	1 ea	Contact for details
116002512	QuickPrep™ 24 x 2 mL Adaptor	1 ea	Contact for details

APPENDIX 1: FASTPREP® INSTRUMENTS

CRYOGENIC ADAPTORS

Our cryogenic Adaptors are ideally suited for extractions of any temperature-unstable or sensitive biological compounds including RNA, siRNA, metabolites, intermediates, and enzymes from even the hardest samples to lyse. Allowing simultaneous cryogenic lysis, the cool adaptors ensure efficient cooling of samples through passive temperature control technology with dry ice. Due to high heat transfer capacity and precise settings of lysis parameters, samples can be repeatedly homogenized without any increase in temperature.

<p>CoolPrep™ 24x2 mL Adaptor Cryogenic adapter holds 24 x 2ml Lysing Matrix tubes.</p>	
<p>CoolTeenPrep™ 6x15 mL Adaptor Cryogenic adapter holds 6 x 15ml Lysing Matrix tubes.</p>	
<p>CoolBigPrep™ 2x50 mL Adaptor Cryogenic adapter holds 2 x 50ml Lysing Matrix tubes.</p>	

Adaptor Ordering Information

Cat No	Item	Unit	Price
116002528	CoolPrep™ 24x2 mL Adaptor	1 ea	Contact for details
116002530	CoolTeenPrep™ 6x15 mL Adaptor	1 ea	Contact for details
116002531	CoolBigPrep™ 2x50 mL Adaptor	1 ea	Contact for details

APPENDIX 1: FASTPREP® INSTRUMENTS

METAL ADAPTORS

Metal Adaptors are ideally suited for work with highly infectious, pathogenic, or other biologically hazardous samples. They withstand temperatures up to 450°C, allowing for sterilization by pyrolysis or autoclaving. Pathogens, including bacteria, viruses, fungi, parasites, viroids, and prions, can thus be effectively eliminated. All-Metal adapters are also safe to use with most laboratory detergents and sterilization solutions, ensuring easy care and maintenance.

<p>QuickPrep™ 24 x 2 mL All Metal Adaptor All-Metal adapter holds 24 x 2ml Lysing Matrix tubes.</p>	
<p>TeenPrep™ 12 x 15 mL All Metal Adaptor Cryogenic adapter holds 12 x 15ml Lysing Matrix tubes.</p>	
<p>BigPrep™ 2 x 50 mL All Metal Adaptor Cryogenic adapter holds 2 x 50 Lysing Matrix tubes.</p>	
<p>Ratchet Nut v3.0 Blue "Locking ratchet highly recommended for all Metal adapter. Alternative locking system for all adapters."</p>	

Adaptor Ordering Information			
Cat No	Item	Unit	Price
116002545	Metal QuickPrep™ 24 x 2 mL All Metal Adaptor	1 ea	Contact for details
116002546	Metal TeenPrep™ 12 x 15 mL All Metal Adaptor	1 ea	Contact for details
116002547	Metal BigPrep™ 2 x 50 mL All Metal Adaptor	1 ea	Contact for details
116005510	Ratchet Nut v3.0 Blue	1 ea	Contact for details

APPENDIX 2: FASTPREP® LYSING MATRIX

APPENDIX 2: FastPrep® Lysing Matrix

Lysing matrices are critical components of the FastPrep® System and help achieve optimal lysing performance. These MP FastPrep® Lysing Matrix Tubes are prepared and dispensed under rigorous conditions, resulting in a premium product that is ready to use with confidence.

Matrices are available as part of MP FastPrep® purification kits, as well as separately. A description of the MP FastPrep® Lysing Matrix Tubes available from MP Biomedicals is included below.

MP FastPrep Lysing Matrix	Description	MP FastPrep Kit (when available)	Sample Type	Extraction Target
A	Contains gamet matrix and ¼-inch ceramic sphere with an orange cap: Use for all sample types except soil.	FastDNA™ SPIN Kit for Plant and Animal Tissue FastDNA™ Kit FastDNA™ Spin Kit	Animal Bacteria Yeast Fungi Plant	DNA RNA Proteins
B	Contains 0.1 mm silica spheres with a blue cap: Use for gram-positive and gram-negative bacteria.	FastRNA™ Pro Blue Kit FastRNA™ SPIN Kit for Microbes FastPROTEIN™ Blue Matrix	Bacteria Spores	DNA RNA Proteins
C	Contains 1-mm silica spheres with a red cap: Use for yeast and fungi.	FastRNA™ Pro Red Kit FastPROTEIN™ Red Matrix	Yeast Fungi	DNA RNA Proteins
D	Contains 1.4 mm ceramic spheres with a green cap: Use for plant and animal tissues.	FastDNA™ SPIN Kit for Plant and Animal Tissue FastRNA™ Pro Green Kit	Animal Plant	DNA RNA Proteins
E	Contains 1.4 mm ceramic spheres, 0.1 mm silica spheres, and one 4 mm glass bead with a purple cap: Use for soil and environmental samples.	FastDNA™ SPIN Kit for Soil FastDNA™ SPIN Kit for Feces FastRNA™ Pro Soil Kits	Soil Sediments Water Feces	DNA RNA Proteins
F	Contains 1.6 mm aluminum oxide particles and 1.6 mm silicon carbide particles with a white cap: Use for breaking tough, hard, or brittle cell membranes from plant, animal, bacteria, mold and other sample types.	FastDNA™ SPIN Kit for Plant and Animal Tissue FastDNA™ Spin Kit FastDNA™ Kit	Plant Animal Bacteria Molds Fungi Coral Emulsions Fixed Samples	DNA RNA Proteins
G	Contains 1.6 mm silicon carbide particles and 2 mm glass beads with a brown cap: Use for breaking tough, hard, or brittle cell membranes from yeast, fungi and spores, as well as brittle plant and animal tissue and more.	FastDNA™ SPIN Kit for Plant and Animal Tissue FastDNA™ Spin Kit FastDNA™ Kit	Plant Animal Yeast Fungi Spores	DNA RNA Proteins
H	Contains 2 mm glass beads and 2 mm yellow zirconium oxide beads with a yellow cap: Use for breaking tough, hard cells including organisms with dense exterior matrices.	FastDNA™ SPIN Kit for Plant and Animal Tissue FastDNA™ Spin Kit FastDNA™ Kit	Plant Animal Wood Seeds Very Dense Soils/Clays	DNA RNA Proteins

APPENDIX 2: FASTPREP® LYSING MATRIX

			Bacterial Aggregates Whole Insects Ticks Ancient /Dried Samples	
I	Contains 2 mm yellow zirconia beads and a 4 mm black ceramic sphere with a clear cap: Use to break primarily by impaction, very tough, hard samples such as chitin exoskeletons and dry grinding of fungal spores	FastDNA™ SPIN Kit for Plant and Animal Tissue FastDNA™ Spin Kit FastDNA™ Kit	Seeds Very dense soils/clays Bacterial Aggregates Whole Insects Ticks Ancient /Dried Samples Dry Grinding Fungal Spores/Rusts	DNA RNA Proteins
J	Contains 2 mm yellow zirconia beads and 1.6 mm aluminum oxide particles and a pink cap: Use for high impaction and low shearing such as isolation of intact organelles and super-molecular structures from tissue, bacteria, molds, and fungi	FastDNA™ SPIN Kit for Plant and Animal Tissue FastDNA™ Spin Kit FastDNA™ Kit	Plant Animal Bacteria (gram +/-) Molds Fungi Coral Emulsions Fixed Samples Fungal Spores /Rusts	DNA RNA Proteins
K	Contains 0.8 mm zirconium silicate beads with a brown cap: Use to break spores, cysts and yeast polysaccharide capsules	GeneClean™ for Ancient DNA Kit	Non-viable Tissues Bone Preserved Tissues Animal By-Products Yeast Spores	DNA RNA Proteins
M	Contains one ¼" (6.35 mm) diameter zirconium oxide coated ceramic grinding sphere, and includes an extra bag of spheres so that 2 can be used to grind very difficult samples by impaction.	FastDNA™ SPIN Kit for Plant and Animal Tissue FastDNA™ Spin Kit FastDNA™ Kit	Tough Tissues Skeletal/Muscle Tissue Lung Tissue Heart Tissue Bone Tissue Seeds Spores	DNA RNA Proteins
S	Contains six (6) 1/8" (3.175 mm) diameter stainless steel beads: Use for RNA and protein extraction from difficult samples where lysis by impaction is preferable.	FastDNA™ SPIN Kit for Plant and Animal Tissue FastDNA™ Spin Kit FastDNA™ Kit	Tough Tissues Skeletal/Muscle Tissue Lung Tissue Heart Tissue Bone Tissue Seeds Spores	DNA RNA Proteins
SS	Contains fifty (50) ¼" (6.35 mm, 0.25 cal) stainless-steel grinding balls: Use for RNA extraction from difficult samples where lysis by impaction is preferable.	FastDNA™ SPIN Kit for Plant and Animal Tissue FastDNA™ Spin Kit FastDNA™ Kit	Tough Tissues Skeletal/Muscle Tissue Lung Tissue Heart Tissue Bone Tissue Seeds Spores	DNA RNA Proteins
Y	Contains 0.5 mm Ytria-Satbilized zirconium oxide spheres with a white cap: Use for yeast, fungi and algal samples.	FastRNA™ SPIN Kit for Yeast	Yeast Fungi Algae	DNA RNA Proteins
Z	Contains 2.0 mm Ytria-Satbilized zirconium oxide spheres with a natural (clear) cap: Use for tough plant and animal samples.	FastDNA™ SPIN Kit for Plant and Animal Tissue	Animal Plant	DNA RNA Proteins

APPENDIX 3: RECOMMENDED PROGRAM SETTINGS

APPENDIX 3: RECOMMENDED PROGRAM SETTINGS

Category: Bacteria							
Preset Name	Speed (m/s)	Adapter	Time (seconds)	Cycles	Rest Time	Lysing Matrix	Quantity (mg)
<i>Listeria monocytogenes</i> cells	6.0	QuickPrep-3 / QuickPrep	30	3	300	B	10 ⁹ cells
<i>Streptococcus pyogenes</i> cells	6.0	QuickPrep-3 / QuickPrep	20	1	N/A	B	10 ⁹ cells
<i>Streptococcus mutans</i> cells	6.0	QuickPrep-3 / QuickPrep	30	1	N/A	B	10 ⁹ cells
<i>Staphylococcus aureus</i> cells	6.0	QuickPrep-3 / QuickPrep	40	2	300	B	10 ⁸ cells
<i>Photobacterium luminescens</i> cells	6.0	QuickPrep-3 / QuickPrep	30	2	300	B	10 ⁹ cells
<i>Escherichia coli</i> cells	6.0	QuickPrep-3 / QuickPrep	30	1	N/A	B	10 ⁸ cells
<i>Mycobacterium tuberculosis</i> cells	6.0	QuickPrep-3 / QuickPrep	45	2	300	B	10 ⁸ cells
<i>Lactococcus lactis</i> cells	6.0	QuickPrep-3 / QuickPrep	30	3	300	B	10 ⁸ cells
Category: Environmental							
Preset Name	Speed (m/s)	Adapter	Time (seconds)	Cycles	Rest Time	Lysing Matrix	Quantity (mg)
Sediment - Soil / Rocks	5.5	QuickPrep-3 / QuickPrep	30	2	300	E	50
Soil - Sandy Sample	4.0	QuickPrep-3 / QuickPrep	30	4	300	E	50
Soil - Litter	5.5	QuickPrep-3 / QuickPrep	30	1	N/A	E	50
Waste Water	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	E	1 mL
Marine Sediment	5.5	QuickPrep-3 / QuickPrep	40	2	300	E	500
Soil from Grassland	5.5	QuickPrep-3 / QuickPrep	30	2	300	E	500

APPENDIX 3: RECOMMENDED PROGRAM SETTINGS

Rhizosphere	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	E	500
Asphalt permeated soil	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	E	500
Category: Feces							
Preset Name	Speed (m/s)	Adapter	Time (seconds)	Cycles	Rest Time	Lysing Matrix	Quantity (mg)
Human Feces	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix E	300
Category: Human / Animal							
Preset Name	Speed (m/s)	Adapter	Time (seconds)	Cycles	Rest Time	Lysing Matrix	Quantity (mg)
Human Lung	6.0	QuickPrep-3 / QuickPrep	30	4	300	Lysing Matrix D	50
Human Breast	6.0	QuickPrep-3 / QuickPrep	30	2	300	Lysing Matrix D	80
Human Kidney	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	50
Human Thyroid Tumors	6.0	QuickPrep-3 / QuickPrep	30	3	300	Lysing Matrix A	100
Mouse Eye	6.0	QuickPrep-3 / QuickPrep	30	4	300	Lysing Matrix D	10
Mouse Heart	6.0	QuickPrep-3 / QuickPrep	30	4	300	Lysing Matrix D	70
Mouse Kidney	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	50
Mouse Femur	6.0	QuickPrep-3 / QuickPrep	30	4	300	Lysing Matrix A	40
Mouse Leg Muscle	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	50
Mouse Intestine	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	50
Mouse Ear	6.0	QuickPrep-3 / QuickPrep	30	4	300	Lysing Matrix D	45
Mouse Tail	6.0	QuickPrep-3 / QuickPrep	30	4	300	Lysing Matrix A	100

APPENDIX 3: RECOMMENDED PROGRAM SETTINGS

Mouse Spleen	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	70
Mouse Lung	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	50
Mouse Liver	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	50
Mouse Brain	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	50
Mouse Pancreatic Cells (bHC9)	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	10 ⁷ cells
Human Ovary Biopsy	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix A	
Human Skin	6.0	QuickPrep-3 / QuickPrep	40	3	300	Lysing Matrix D	19
Category: Plant							
Preset Name	Speed (m/s)	Adapter	Time (seconds)	Cycles	Rest Time	Lysing Matrix	Quantity (mg)
Alpowa Wheat Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	75
Alpowa Wheat Seed	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix A	100
<i>Arabidopsis thaliana</i> Fresh Leaves (50 mg)	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	50
<i>Arabidopsis thaliana</i> Fresh Leaves (200 mg)	6.0	QuickPrep-3 / QuickPrep	40	2	300	Lysing Matrix D	200
Bartlett Pear Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	50
Classic Oat Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	75
Classic Oat Seed	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix A	100
Corn Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	100
Crest Barley Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	100
Crest Barley Root	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix A	300

APPENDIX 3: RECOMMENDED PROGRAM SETTINGS

Kaybonnet Rice Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	100
Kaybonnet Rice Seed	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix A	100
Klages Barley Root	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix A	300
Klages barley Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	75
Tobacco Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	75
Lafitte Rice Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	2	300	Lysing Matrix D	75
Lafitte Rice Sprout Leaf	6.0	QuickPrep-3 / QuickPrep	30	1	N/A	Lysing Matrix D	100
Soybean Seed	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix A	100
Corn Seed	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix A	100
Oat FL 502 Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	75
Oact FL 502 Seed	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix A	100
Riser Oat Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	70
Richland Soybean Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	100
Tam Wheat Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	75
Tam Wheat Root	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix A	80
Tomato, Early Girl, Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	30	4	300	Lysing Matrix D	75
Williams 82 Soybean Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	70
Wrens Rye Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	100
Pine Needle	6.0	QuickPrep-3 / QuickPrep	30	1	N/A	Lysing Matrix A	100
Basil Leaf Tissue	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix D	50
Cassava Root	6.0	QuickPrep-3 / QuickPrep	60	1	N/A	Lysing Matrix A	80

APPENDIX 3: RECOMMENDED PROGRAM SETTINGS

Category: Yeast & Fungi							
Preset Name	Speed (m/s)	Adapter	Time (seconds)	Cycles	Rest Time	Lysing Matrix	Quantity (mg)
<i>Saccharomyces cerevisiae</i> cells	6.0	QuickPrep-3 / QuickPrep	40	1	N/A	Lysing Matrix	10 ⁸ cells
<i>Schizosaccharomyces pombe</i> cells	5.0	QuickPrep-3 / QuickPrep	15	4	300	Lysing Matrix	10 ⁸ cells
<i>Candida albicans</i> cells	6.0	QuickPrep-3 / QuickPrep	30	2	300	Lysing Matrix	10 ⁸ cells
<i>Cryptococcus neoformans</i> cells	6.0	QuickPrep-3 / QuickPrep	30	4	300	Lysing Matrix	10 ⁸ cells
<i>Aspergillus fumigatus</i> cells	6.0	QuickPrep-3 / QuickPrep	30	2	300	Lysing Matrix	10 ⁸ cells
<i>Fusarium Solani</i> cells	6.0	QuickPrep-3 / QuickPrep	30	2	300	Lysing Matrix	10 ⁸ cells

APPENDIX 4: FASTPREP® INSTRUMENTS FOR OTHER APPLICATIONS

FastPrep-96™ Instrument**Nothing Resists Grinding, Lysis and Homogenization with the High-Throughput FastPrep-96™ System**

MP Biomedicals' FastPrep-96™ ([Figure A1](#)) is a versatile high-speed homogenizer offering the ultimate in high-throughput sample preparation. Developed for resistant samples, the FastPrep-96™ instrument uses a high-speed linear motion to disrupt thoroughly any tissues and cells in just seconds through the simultaneous beating of specialized Lysing Matrix beads on the sample material. This high-performance system saves hours of work during the sample preparation stage.

- High Throughput :** Process up to 192 samples simultaneously in 2 x 96 deep well plates
- Excellent Reproducibility:** Consistent lysis results in every well with automated feedback control
- Fast Processing Speed:** 1800 Oscillations/min and 1.5-inch stroke — highest available
- True Linear Motion:** Eliminates the need to re-orient plates mid-cycle



2 x 96 well plate adapter

included with instrument



Figure A1: FastPrep-96™ Instrument

The FastPrep-96™ instrument offer the highest power settings and widest variety of Adaptors available (2 x 96 deep well plates, 96 x 2 mL, 48 x 4.5 mL, 24 x 15 mL, 8 x 50 mL, and 1 x 250 mL tubes) ([Figure A2](#)) and with the closed-loop, automated control of lysing power and time, the new FastPrep-96™ is the perfect solution for all of your sample preparation needs.

APPENDIX 4: FASTPREP® INSTRUMENTS



Figure A2: FastPrep-96 Adaptors

FastPrep-96™ Ordering Information			
Cat No.	Item	Unit	Price
116010500	FastPrep-96™ Instrument (includes 2 x 96 well plate adapter)	1 ea	Contact for details
116010570	QuickFlex™ Adaptor (96 x 2 mL)	1 ea	Contact for details
116010580	TallFlex™ Adaptor (48 x 4.5 mL)	1 ea	Contact for details
116010560	TeenFlex™ Adaptor (24 x 15 mL)	1 ea	Contact for details
116010550	BigFlex™ Adaptor (8 x 50 mL)	1 ea	Contact for details
116010590	LargeFlex™ Adaptor (2 x 250 mL)	1 ea	Contact for details
116010595	ConeFlex™ Legacy Adapter (allows any existing FastPrep-24™ adapters to be used on the FastPrep-96™ instrument)	1 ea	Contact for details

APPENDIX 4: FASTPREP® INSTRUMENTS

SuperFastPrep-2™ Portable Homogenizer, 2 x 2 mL, battery operated

The Fastest and Most Powerful Handheld SamplePrep Tool

- Omnidirectional motion with the highest G's!
- Handheld, cordless sample preparation instrument for lab and field use.
- Uses standard 2 mL screw cap tubes for lysis.
- Compatible with all FastPrep® Lysing Matrices.



Figure A3: The SuperFastPrep-2™

The SuperFastPrep-2™ ([Figure A3](#)) from MP Biomedicals is the newest addition to the trusted FastPrep Instruments for sample grinding. For the first time, FastPrep® technology is available in a lightweight, compact, handheld format. An innovation in the sample lysis industry, the SuperFastPrep-2™ is a portable omnidirectional bead beating system that attaches to a rechargeable, battery-powered rotary tool.

SuperFastPrep-2™ Ordering Information			
Cat No.	Item	Unit	Price
116012500	SuperFastPrep-2™ Instrument	1 ea	Contact for details

Worldwide Headquarters

MP Biomedicals, LLC
9 Goddard
Irvine, CA
92618
USA
Contact number: 1.800.854.0530
custserv.na@mpbio.com

North America

MP Biomedicals North America

MP Biomedicals
NA Distribution & Customer Service Center
29525 Fountain Parkway
Solon, OH
44139
USA
Contact Number: +1 800.854.0530
Fax No: +1 440.337.1180
custserv.na@mpbio.com

South America

MP Biomedicals Brasil

MP Biomedicals do Brazil LTDA
Rua Cavalheiro Ernesto Giuliano, 909
São Caetano do Sul- SP
CEP 09540-400
BRAZIL
Contact Number: +55 11 3565-7187
custserv.la@mpbio.com

Europe, Middle East, Africa

MP Biomedicals France

MP Biomedicals SARL
Parc d'Innovation BP50067
Rue Geiler de Kaysersberg
Illkirch Cedex
67402
FRANCE
Contact Number: 00800.7777.9999 or +33 3.88.67.19.45
Fax No: +33 3.88.67.19.45
custserv.fr@mpbio.com

MP Biomedicals Germany

MP Biomedicals Germany GmbH
Thüringer Straße 15
Eschwege
37269
GERMANY
Contact Number: 0800.426.67.337 or 00800.7777.9999
Fax No: 0800.629.67.337
custserv.de@mpbio.com

MP Biomedicals Serbia

MP Bio d.o.o.
Prokupacka 41
Belgrade
11000
SERBIA
Contact Number: +381 11 2648 236 or +381 11 2648 148
Fax No: +381 11.242.1949
custserv.se@mpbio.com

MP Biomedicals Russian Federation

MPBA Diagnostics LLC (MP Biomedicals Russia)
3 Marksistskaya St.
Bldng 2
Moscow
109147
RUSSIAN FEDERATION
Contact Number: +7 495 604-13-44
Fax No: +7 495 604-13-44
custserv.rs@mpbio.com

Asia

MP Biomedicals Asia Pacific

MP Biomedicals Asia Pacific Pte Ltd, HQ
2 Pioneer Place
627885
SINGAPORE
Contact Number: +65 6775.0008
Fax No: +65 6774.6146
custserv.ap@mpbio.com

MP Biomedicals Japan

MP Bio Japan K. K.
Seven Bldg. 4F, 3-7-14
Nihonbashi-Ningyocho
Chuo-ku, Tokyo
103-0013
JAPAN
Contact Number: +81 3.6667.0730
Fax No: +81 3.6667.0733
custserv.jp@mpbio.com

MP Biomedicals China

MP Biomedicals (Shanghai) Co., Ltd.
Unit 02, 2nd floor, Building #6
1000 Jinhai Road
Pudong District
Shanghai
People's Republic of CHINA
Contact Number: +86 400.1500.680
Fax No: +86 21.6258.6278
custserv.cn@mpbio.com

MP Biomedicals India

MP Biomedicals (India) Pvt. Ltd
Plot No C -57/3, TTC Industrial Area
Pawane Village, Turbhe, Navi Mumbai - 400 705
Mumbai
400705
INDIA
Contact Number: +91 22.27636921/22/25
custserv.in@mpbio.com

MP Biomedicals Korea

MP Biomedicals Korea
B-809 Garden Five Works,
52, Chungmin-ro,
Songpa-gu, Seoul,
05839
Korea
Contact Number: +82 2.425.5991
Fax No: +82 2.425.5997
custserv.kr@mpbio.com

Australia, New Zealand

MP Biomedicals Australia

MP Biomedicals Australia Pty Ltd
P.O. Box 187
2 / 29 Bearing Road
Seven Hills
NSW 2147
Australia
Contact Number: 1800249998 or +61 2.8824.2100
Fax No: +61 2.9838.7390
custserv.au@mpbio.com

MP Biomedicals New Zealand

MP Biomedicals New Zealand
37-39 Waipareira Avenue
Henderson
Auckland
0610
New Zealand
Postal address:
P.O. Box 1607
Shortland Street
Auckland
1140
New Zealand
Contact Number: +64 9.912.246
Fax No: +64 9.838.4209
custserv.nz@mpbio.com