

SPINeasy RNA Kit for Bacteria (with Lysing Matrix)

For the isolation of RNA from bacteria samples

Size: 50 & 5 preps

Storage: 15-25 °C

Cat. No.: 116541050 (50 PREPS)

116541000 (5 PREPS)

Content Version: November 2023



Table of Contents

1. Introduction to SPINeasy RNA Kit for Bacteria	3
2. Kit Components and User Supplied Materials	4
3. Storage and Kit Stability	5
4. Important Consideration Before Use	5
5. Safety Precautions	6
6. Protocol.....	7
7. Flow Chart.....	10
8. Data	11
9. Troubleshooting	13
10. Product Use Limitation & Warranty	14

1. Introduction to SPINeasy RNA Kit for Bacteria

SPINeasy RNA Kit for Bacteria is a silica-membrane spin-column kit that enables quick and convenient purification of total RNA from gram-positive and gram-negative bacteria. Included in the kit is our specially formulated RNASS solution that stabilizes and protects RNA in bacteria samples. The use of Lysis Buffer R and Lysing Matrix B in combination with FastPrep® instruments from MP Biomedicals enables highly efficient lysis of bacterial samples within seconds. With a simple workflow, the kit allows multiple samples to be processed simultaneously, without the use of toxic substances such as phenol and chloroform. Total RNA of high quality and integrity can be purified typically within 40 minutes and is immediately available for RT-PCR and other downstream applications.

Visit www.mpbio.com to explore additional products to support your research.

Kit Specifications at a Glance

Technology	Silica membrane technology
Format	Spin column
Sample	Bacteria (e.g., gram-positive/negative bacteria)
Sample amount	up to 6 mL bacterial culture
Typical yield	up to 800 ng/μL (sample dependent)
Elution volume	50-100 μL
Preparation time	<40 min

2. Kit Components and User Supplied Materials

2.1 SPINeasy RNA Kit for Bacteria Component

Components	50 PREPS (Cat.No.: 116541050)		5 PREPS (Cat.No.: 116541000)	
	Package	Cat. No.	Package	Cat. No.
Equilibration buffer R	12 mL	112075704	1.2 mL	112075804
Lysing Matrix B	50 ea	116911050	5 ea	116911005
RNASS	15 mL	116541051	1.5 mL	116541001
Lysis Buffer R	60 mL	116541052	6.0 mL	116541002
Wash Buffer R	6 mL	116541053	600 µL	116541003
Nuclease-free water	10 mL	116541054	100 µL	116541004
DNase I	1 vial	116541055	1 vial	116541055
DNase I Buffer	5 mL	116541056	500 µL	116541006
Column R with collection tube	50 ea	116541057	5 ea	116541007
Gram-positive Bacteria Pre-treatment Buffer	12 mL	116533051	1.2 mL	116533001
Quick-start protocol	1 ea	-	1 ea	-
Instruction Manual	Available www.mpbio.com			
MSDS & CoA	Available www.mpbio.com			

2.2 User Supplied Materials

- FastPrep® Instrument - FastPrep-24™ 5G (Cat. No.116005500) or Vortex.
- Microcentrifuge capable of spinning at $\geq 14,000 \times g$.
- Absolute ethanol (50 mL for preparing Wash Buffer R and 750 µL per prep for sample preparation)
- Nuclease-free 2 mL microcentrifuge tubes
- Nuclease-free 1.5 mL microcentrifuge tubes
- Optional: To prepare 20 mg/mL lysozyme (to be supplied by user) with Gram-positive Bacteria Pre-treatment Buffer included in the kit. Store at 4 °C after adding lysozyme.

3. Storage and Kit Stability

DNase I should be stored at 2-8 °C. All other components and reagents of the SPINeasy RNA Kit for Bacteria can be stored at room temperature (15-25 °C) until the expiration date printed on the kit label. For extended storage or storage in dry condition (humidity < 40%), store the columns at 2-8 °C to maintain performance.

4. Important Consideration Before Use

- Add 50 mL (5 mL for sample kit) of absolute ethanol to Wash Buffer R and mark the bottle.
- Prepare DNase I solution according to instructions in Section 6.1.
- Prepare one 2 mL microcentrifuge tube per prep for lysate binding preparation.
- Prepare one 1.5 mL microcentrifuge tube per prep for elution of purified RNA.
- Prepare one 1.5 mL microcentrifuge tube for DNase I-buffer mixture.
- Lysis can be performed by vortexing the sample in a vial of Lysing Matrix B at the maximum speed if a FastPrep instrument is unavailable.
- Centrifugation speed stated in the manual will be a guideline; use the maximum speed available if 14,000 x g is not feasible.

5. Safety Precautions

Lysis Buffer R contains a component that can be harmful if swallowed and may cause irritation when in contact with skin and eyes. To prevent accidental ingestion, do not eat, drink or smoke when using this product. Wear personal protective equipment (gloves, lab coat and eye protection) to prevent contact with the skin or mucous membranes. Consult the Material Safety Data Sheet at www.mpbio.com for additional details.

6. Protocol

Column Preparation

Pipette 200 μL Equilibration Buffer into Column MB. Incubate for 1 min at room temperature and centrifuge for 30 sec @ maximum speed. Discard flow-through and reuse the collection tube.

6.1 Preparation of DNase I solution

Briefly spin down the vial of lyophilized DNase I provided and resuspend with 500 μL of Nuclease-free water. Mix well to dissolve. Store DNase I solution at $-20\text{ }^{\circ}\text{C}$ in aliquots and avoid repeated freeze-thawing. Note: Do not prepare DNase I solution in DNase I Buffer.

6.2 Pre-treatment of Gram-positive Bacteria (Optional)

If necessary, pre-treatment of gram-positive bacteria with lysozyme may be performed, which may result in higher RNA yields. Good quality RNA of reasonable yields can also be obtained without performing the pre-treatment.

1. To prepare 20 mg/mL lysozyme (to be supplied by user) with Gram-positive Bacteria Pre-treatment Buffer included in the kit. Store at $4\text{ }^{\circ}\text{C}$ after adding lysozyme.
2. Harvest cells by centrifugation at $10,000 \times g$ for 3 mins and discard supernatant. Resuspend bacterial pellet in 200 μL of lysozyme pre-treatment buffer.
3. Incubate for 30 mins at $37\text{ }^{\circ}\text{C}$.
4. Transfer sample to a vial of Lysing Matrix B, add 250 μL of RNASS and 750 μL of Lysis Buffer R.
5. Proceed to step 6.3 of the RNA Extraction Protocol.
6. Note that in this case, up to 950 μL of the lysate supernatant may be used in step 6 of the RNA Extraction Protocol. Consequently, add 1:1 volume of ethanol proportionately and repeat sample loading steps (steps 8 and 9) to load all the mixture.


6.3 RNA Extraction Protocol

1. Pellet 1 - 6 mL bacterial culture by centrifugation at $10,000 \times g$ for 3 mins and discard supernatant
2. Resuspend cell pellet in 250 μL of RNASS and transfer to a vial of Lysing Matrix B.

3. Add 750 μ L of Lysis Buffer R.
4. Homogenize in a FastPrep instrument for 40 seconds at speed setting of 6.0 m/s. Alternatively, vortex samples at the maximum speed for 5 mins if a FastPrep instrument is not available. It is recommended to perform vortexing with the use of an adapter to hold the vials, especially if multiple samples are to be processed simultaneously.
5. Centrifuge at 14,000 x g for 10 mins. Note: Centrifuge at the maximum speed for all steps if 14,000 x g is not feasible.
6. Carefully pipette out 750 μ L of the supernatant to a clean, nuclease-free 2 mL microcentrifuge tube.
7. Add 1:1 volume of absolute ethanol (e.g., 750 μ L ethanol to 750 μ L lysate supernatant) and mix well by pipetting up and down.
8. Transfer 750 μ L of the mixture to a Column R with collection tube.
9. Centrifuge at 14,000 x g for 1 min. Discard flow through and reuse collection tube.
10. Repeat steps 8 and 9 to load the remaining mixture.
11. DNase I digestion: In a clean 1.5 mL microcentrifuge tube, add 5 μ L of DNase I solution to 75 μ L of DNase I Buffer per prep. Mix well and add 80 μ L to the center of the column membrane. Incubate at room temperature for 15 mins. Note: Mixture of DNase I solution with DNase I buffer should be freshly prepared prior to usage. It is not recommended to store the remaining mixture.
12. Add 500 μ L of Wash Buffer R to the column.
13. Centrifuge at 14,000 x g for 1 min. Discard flow through and reuse collection tube.
14. Add 500 μ L of Wash Buffer R to the column.
15. Centrifuge at 14,000 x g for 1 min. Discard flow through and reuse collection tube.
16. Centrifuge at 14,000 x g for an additional 1 min to dry column.
17. Remove collection tube and place column onto a clean 1.5 mL microcentrifuge tube.
18. Add 100 μ L of Nuclease-free water to the center of the membrane. Incubate at room temperature for 1 min. For samples with low RNA content, reducing the elution volume to 50 μ L may increase the concentration of eluted RNA.

19. Centrifuge at 8,000 x g for 1 - 2 mins to elute RNA.
20. Eluted RNA will be collected in the microcentrifuge tube. For best results, proceed to perform downstream applications immediately and keep RNA chilled on ice while working to prevent degradation. Store remaining RNA at -80 °C in aliquots and avoid repeated freeze-thawing.

7. Flow Chart

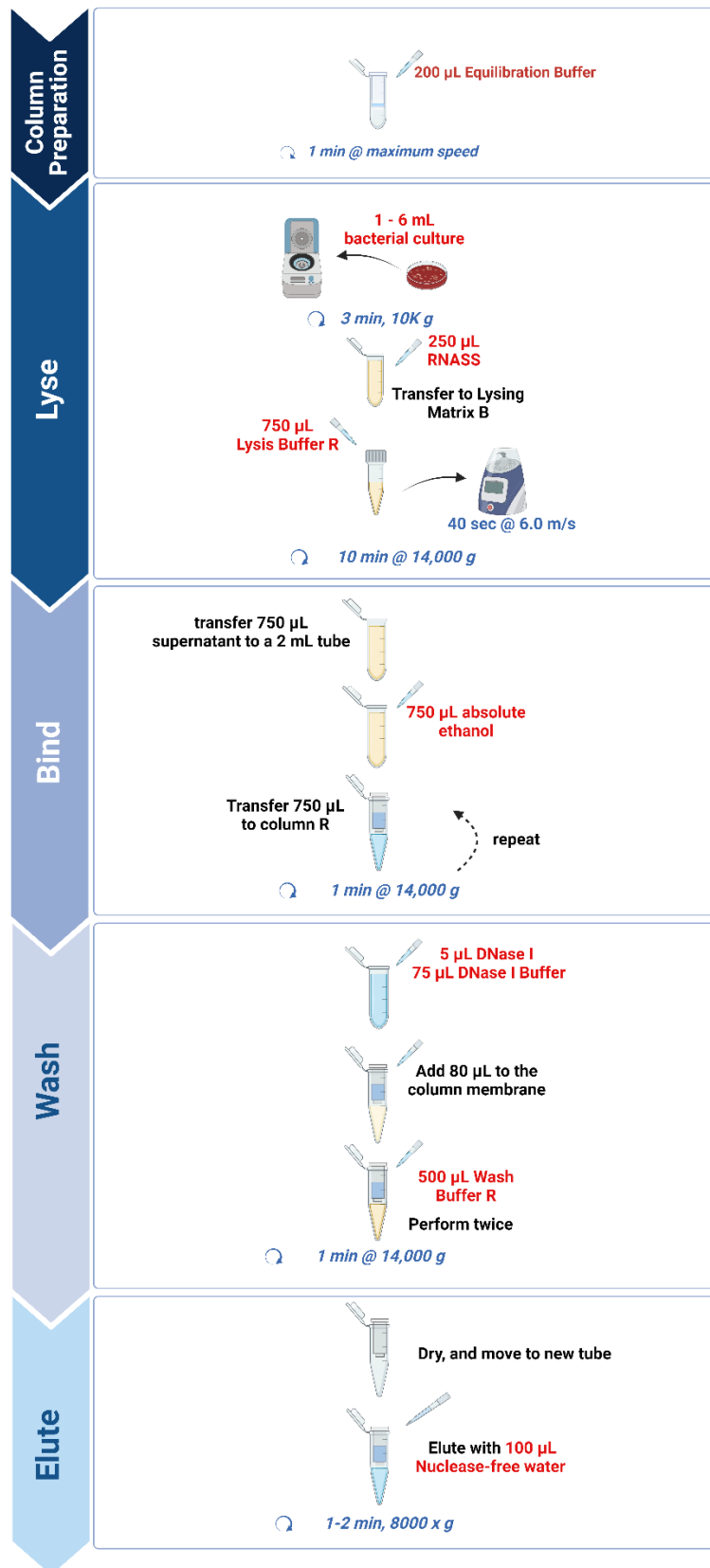

8. Data

The following results demonstrate examples of RNA extracted from bacteria samples using SPINeasy RNA Kit for Bacteria. Extracted RNA are of high yield, quality and integrity, as indicated by concentration readings, A₂₆₀/A₂₈₀ and A₂₆₀/A₂₃₀ ratios of around 2, and 23S and 16S bands clearly resolved in agarose gel electrophoresis. RNA extracted with this kit can be amplified by RT-PCR.

Table 1. Quality and quantity of RNA extracted from representative bacteria samples using SPINeasy RNA Kit for Bacteria.

Sample	Starting Amount	Extraction Results		
		Concentration	A ₂₆₀ /A ₂₈₀	A ₂₆₀ /A ₂₃₀
<i>E. Coli</i>	Pellet from 3 mL overnight culture	763.59	2.10	2.26
<i>S. aureus</i>	Pellet from 3 mL overnight culture	349.47	2.13	2.04

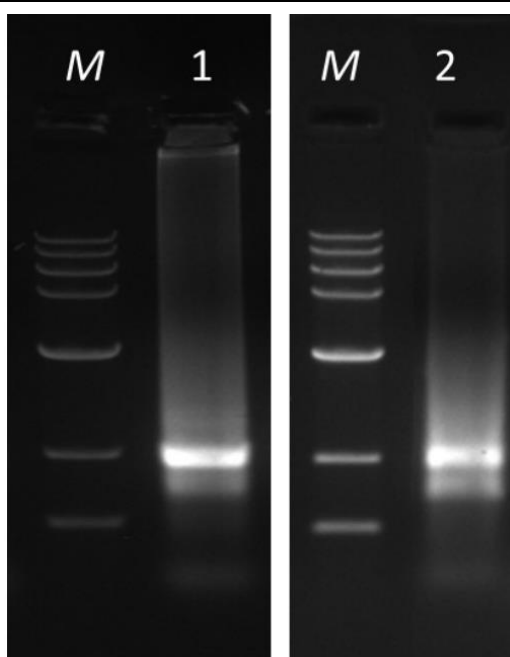


Figure 1: Total RNA extracted from bacteria samples using SPINeasy RNA Kit for Bacteria, analyzed using agarose gel electrophoresis. M: DNA marker; Lane 1: *E. coli*; Lane 2: *S. aureus*.

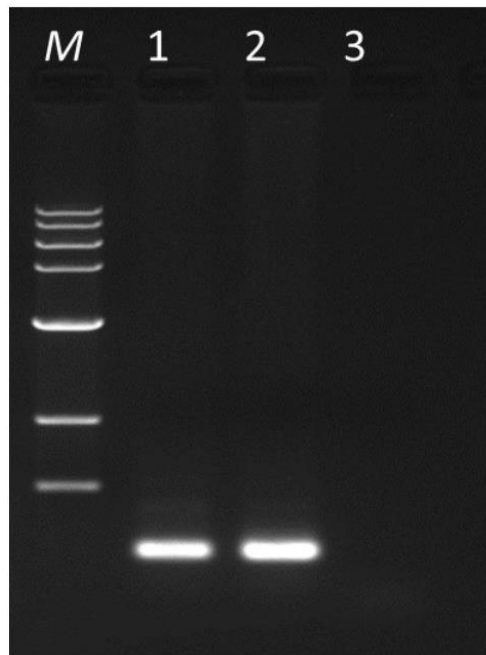


Figure 2: RT-PCR amplification of RNA extracted from bacteria samples using SPINeasy RNA Kit for Bacteria. M: DNA marker; Lane 1: *E. coli*; Lane 2: *S. aureus*; Lane 3: Negative control.

9. Troubleshooting

This guide may be useful in solving any problems that may arise. For further assistance, please contact our technical support team at apac-techsupport@mpbio.com

Problem	Recommendation
Low RNA Yield / reduced RNA integrity	Ensure the extraction was performed according to kit manual instructions.
	Sample with low RNA content: (i) Increase amount of starting material; (ii) Elute in a smaller volume (50 μ L).
	Insufficient lysis or over-lysis: When using a vortex instead of a FastPrep for sample homogenization, lysis condition may be optimized by testing reduced or extended vortexing duration.
	For gram-positive bacteria, performing pre-treatment with lysozyme can lead to increased RNA yields (Refer to 6.2).
	Sample incompatible with RNASS. While RNA yield and quality are greatly enhanced with the use of RNASS for most bacteria samples, RNASS may not be suitable for use with certain sample types. In these cases, resuspend bacterial pellet with 1 mL of Lysis Buffer R, transfer to Lysing Matrix B and proceed to step 4 of the RNA Extraction Protocol without using RNASS.
	Poor sample quality. For best results, freshly prepared samples should be used.
Smear RNA Bands	RNase contamination. Work with nuclease-free tubes and pipette tips. Handle samples and perform all steps with clean gloves. Decontaminate work surfaces with RNase Erase® (Cat. No. 112440204).
	Poor sample quality. For best results, freshly prepared samples should be used.
	Sample over-lysis. Reduce FastPrep speed and/or duration.
DNA Contamination	RNA degradation. Work with freshly purified RNA and keep RNA chilled on ice after elution. RNA should be stored at -80 °C and freeze-thawing should be avoided.
	Perform on-column DNase I digestion according to step 11 of the RNA extraction protocol. Ensure that DNase I solution is prepared according to manual instructions. Once dissolved, DNase I should be stored at -20 °C.

10. Product Use Limitation & Warranty

The products presented in this instruction manual are for research or manufacturing use only. They are not to be used as drugs or medical devices to diagnose, cure, mitigate, treat, or prevent diseases in humans or animals, either as part of an accepted course of therapy or in experimental clinical investigation. These products are not to be used as food, food additives or general household items. Purchase of MP Biomedicals products does not grant rights to reproduce, modify, or repackage the products or any derivative thereof to third parties. MP Biomedicals makes no warranty of any kind, expressed or implied, including merchantability or fitness for any particular purpose, except that the products sold will meet our specifications at the time of delivery.

Buyer's exclusive remedy and the sole liability of MP Biomedicals hereunder shall be limited to, at our discretion, no replacement or compensation, product credits, refund of the purchase price of, or the replacement of materials that do not meet our specification. By acceptance of the product, Buyer indemnifies and holds MP Biomedicals harmless against, and assumes all liability for, the consequence of its use or misuse by the Buyer, its employees, or others, including, but not limited to, the cost of handling. Said refund or replacement is conditioned on Buyer notifying within thirty (30) days of receipt of product. Failure of Buyer to give said notice within thirty (30) days shall constitute a waiver by the Buyer of all claims hereunder with respect to said material(s).

Australia

Tel: +61 2.8824.2100
Tel: +61 1800.249.998
Email: custserv.au@mpbio.com

Austria & Germany

Tel: 0800.426.67.337
Tel: 00800.7777.9999
Email: custserv.de@mpbio.com

Belgium

Tel: 00800.7777.9999
Email: custserv.be@mpbio.com

Canada

Tel: +1 800.854.0530
Email: custserv.ca@mpbio.com

China

Tel: +86 400.150.0680
Email: custserv.cn@mpbio.com

Europe

Tel: +33 3.88.67.54.25
Tel: +33 00800.7777.9999
Email: custserv.eur@mpbio.com

France

Tel: +33 3.88.67.54.25
Email: custserv.fr@mpbio.com

India

Tel: +91 22.27636921/22/25
Email: custserv.in@mpbio.com

Italy

Tel: 00800.7777.9999
Email: custserv.it@mpbio.com

Japan

Tel: +81 3.6667.0730
Email: custserv.jp@mpbio.com

Latin America

Tel: +1 800.854.0530
Tel: +1 440.337.1200
Email: custserv.la@mpbio.com

New Zealand

Tel: +64 9.912.2460
Email: custserv.nz@mpbio.com

North America

Tel: +1 800.854.0530
Tel: +1 440.337.1200
Email: custserv.na@mpbio.com

Poland

Tel: 00800.7777.9999
Email: custserv.po@mpbio.com

Russia

Tel: +7 495 604.13.44
Email: custserv.rs@mpbio.com

Serbia

Tel: +381 11.242.1972
Email: custserv.se@mpbio.com

Singapore/ APAC

Tel: +65 6775.0008
Tel: +65 6394.7675
Email: custserv.ap@mpbio.com

South Korea

Tel: +82 2.425.5991
Email: custserv.kr@mpbio.com

Switzerland

Tel: 00800.7777.9999
Email: custserv.ch@mpbio.com

The Netherlands

Tel: 00800.7777.9999
Email: custserv.nl@mpbio.com

United Kingdom

Tel: 0800.282.474
Email: custserv.uk@mpbio.com

www.mpbio.com

