

SPINeasy™ Host Depletion Microbial DNA Kit

Cat. No.: 116545050 (50 PREPS) / 116545000 (5 PREPS)



Scan QR code for more information
from instruction manual

Quick-Start Protocol

Revision 2.0 Aug 2023

Notes before starting:

- Upon receipt, store **Host Depletion Enzyme HDE** and **Microbial Selection Buffer** at 4 °C.
- Add 50 mL (5 mL for sample kit) of absolute ethanol to Wash Buffer HD2 and mark the bottle.
- Prepare Host Depletion Enzyme HDE solution: Spin down components briefly and add 120 µL of **HDE Reconstitution Buffer** to the entire content in the vial of **Host Depletion Enzyme HDE**. Mix well to dissolve and store reconstituted solution at -20 °C.

Host Cell Lysis

- 1 In a 2 mL microcentrifuge tube, add 200 µL of sample to 1 mL of **Host Lysis Buffer**.
- 2 Vortex at for 10 seconds.
- 3 Incubate for 15 mins at room temperature with constant end-over-end tube rotation.
- 4 Vortex at for 10 seconds and centrifuge at 10,000 x g for 5 mins.
- 5 Carefully aspirate and discard supernatant without disturbing the pellet.

Host DNA Depletion

- 6 Resuspend pellet in 100 µL of **Host Depletion Buffer**.
- 7 Add 1 µL of Host Depletion Enzyme HDE solution* and vortex briefly to mix.
*Prepare Host Depletion Enzyme HDE solution according to instructions in "Notes before starting"
- 8 Incubate at 37 °C for 10 mins.
- 9 Add 200 µL of **Microbial Selection Buffer** and vortex briefly to mix.
- 10 Incubate at 37 °C for 30 mins.
- 11 Vortex for 10 seconds and spin down briefly to collect contents at the bottom of the tube.

Microbial Cell Lysis

- 12 Add 700 µL of **Microbial Lysis Buffer**.
- 13 Mix by pipetting up and down several times and transfer all the mixture to a vial of **Lysing Matrix E**.
- 14 Homogenize in a FastPrep Instrument for 10 seconds at speed setting of 6.0 m/s.
If a FastPrep Instrument is not available, vortex samples at the maximum speed for 1 – 5 mins.
- 15 Centrifuge at 14,000 x g for 10 mins.

Microbial DNA Purification

- 16 Add 200 µL of **Equilibration Buffer** to the **Column HD with collection tube** membrane. Wait for 1 min and centrifuge at 14,000 x g for 1 min, discard flow through and reuse collection tube.
- 17 Transfer 750 µL of the supernatant to a Column HD with collection tube.
- 18 Centrifuge at 14,000 x g for 1 min. Discard flow through and reuse collection tube.
- 19 Add 500 µL of **Wash Buffer HD1**. Centrifuge at 14,000 x g for 1 min, discard flow through and reuse collection tube.
- 20 Add 750 µL of **Wash Buffer HD2**. Centrifuge at 14,000 x g for 1 min, discard flow through and reuse collection tube.
- 21 Centrifuge at 14,000 x g for an additional 1 min to dry column.
- 22 Remove collection tube and place column onto a clean 1.5 mL microcentrifuge tube.
- 23 Add 50 µL of **Elution Buffer HD** to the center of the membrane. Incubate at room temperature for 5 mins.
- 24 Centrifuge at 8,000 x g for 2 mins to elute microbial DNA. Store DNA at -20 °C.




MP Biomedicals nucleic acid extraction kits are designed for simple, efficient, and rapid purification of DNA and RNA from various types of samples. Our wide range of instruments and reagent kits provide you a one-stop solution for your sample preparation works.



FastPrep® Homogenizer
Powerful, Flexible, Intuitive



Lysing Matrix
Complete, Quantitative lysing method

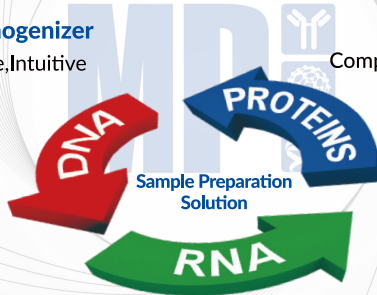

MagBeads FastDNA / FastRNA Kits & SPINeasy™ Extraction Kits
Highly purified DNA / RNA / Protein



Automated Nucleic Acid Purification (aNAP) Instruments
MPure-96™ Process up to 96 samples
MPure-32™ Process up to 32 samples

